


# **Molecular Imager<sup>®</sup> Gel Doc<sup>™</sup> XR+ and ChemiDoc<sup>™</sup> XRS+ Systems with Image Lab<sup>™</sup> Software**

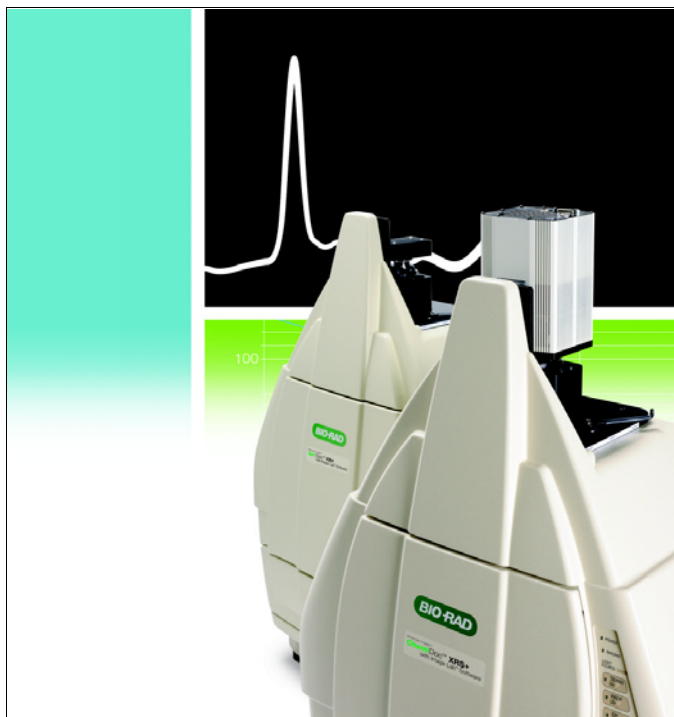



Copyright ©2010 Bio-Rad Laboratories, Inc. Reproduction in any form, either print or electronic, is prohibited without written permission of Bio-Rad Laboratories, Inc.

## **CREDITS**

1. Image Lab software is based in part on the work of the Qwt project (<http://qwt.sf.net>).
2. Image Lab software is based in part on the work of the CImg project (<http://cimg.sourceforge.net/>).  
See license for detail at  
[http://www.cecill.info/licences/Licence\\_CeCILL-C\\_V1-en.html](http://www.cecill.info/licences/Licence_CeCILL-C_V1-en.html)
3. Image Lab software is based in part on the work of the Independent JPEG Group (<http://www.ijg.org/>)

## **TRADEMARKS**

Molecular Imager®, Gel Doc™ XR+, ChemiDoc™ XRS+, and Image Lab™ are trademarks belonging to Bio-Rad Laboratories, Inc.

Windows, Windows XP, Windows Vista, Windows 7, Excel, Microsoft, and Power Point are trademarks of Microsoft Corporation. FireWire, Mac, and Mac OS X are trademarks of Apple Inc. Intel Core is a trademark of Intel Corporation. Hoechst is a trademark of Hoechst Aktiengesellschaft, Mitsubishi is a trademark of Mitsubishi Corporation. SYBR Safe, SYBR Green, SYBR Gold, and SYPRO are trademarks of Invitrogen Corporation. Coomassie is a trademark of BASF Aktiengesellschaft. Cy is a trademark of GE Healthcare group companies. Fluor and Quasar are trademarks of Biosearch Technologies, Inc. Other brands or product names are trademarks of their respective holders.

## **LEGAL NOTICES**

Bio-Rad Laboratories, Inc. is licensed by Invitrogen Corporation to sell SYPRO products for research use only, under US patent 5,616,502.

IL v. 3.0



# Safety and Regulatory Compliance

---

## Important Safety Information

Please read these instructions before attempting to operate the Molecular Imager® Gel Doc™ XR+ or ChemiDoc™ XRS+ imaging systems.

This instrument is suitable for research use only. It must be used, therefore, only by specialized personnel who know the health risks associated with the reagents that are normally used with this instrument.

Use of the Gel Doc XR+ and ChemiDoc XRS+ systems involves UV illumination. Proper precautions must be taken to avoid eye and skin exposure to the UV radiation. This instrument is meant for use only by trained personnel who know the health risks associated with UV radiation normally used with this instrument. The acrylic shield provides some UV protection. However, it does not guarantee complete protection, and it is designed to shield only the person working in front of the system.

**WARNING!** Use of the acrylic screen does not guarantee the user protection from UV radiation. The use of protective eyeglasses, mask, and/or gloves is strongly recommended.

## Warranty

The Gel Doc XR+ and ChemiDoc XRS+ imaging systems are warranted against defects in materials and workmanship for 1 year. If any defect occurs in the instrument during this warranty period, Bio-Rad Laboratories, Inc. will repair or replace the defective parts at its discretion without charge. The following defects, however, are specifically excluded:

- Defects caused by improper operation
- Repair or modification done by anyone other than Bio-Rad Laboratories, Inc. or their authorized agent
- Use of spare parts supplied by anyone other than Bio-Rad Laboratories, Inc.
- Damage caused by accident or misuse
- Damage caused by disaster
- Corrosion caused by improper solvents or samples



## General Precautions

- Please read the instruction manual carefully
- The instrument must be used only for the intended purpose of gel documentation in research laboratories
- The instrument must be connected to a grounded power source line and protected by a circuit breaker
- Do not pour liquids directly on or inside the instrument
- Switch off all the lights immediately after use
- Clean the transilluminator platen after use

## Regulatory Notices

The Gel Doc XR+ and ChemiDoc XRS+ imaging systems are designed and certified to meet EN61010, the internationally accepted electrical safety standards, EMC regulations, and TUV requirements. Certified products are safe to use when operated in accordance with this system instruction manual. This instrument should not be modified or altered in any way. Modification or alteration of this instrument will:

- Void the manufacturer's warranty
- Void the regulatory certifications
- Create a potential safety hazard

**WARNING!** Bio-Rad Laboratories, Inc. is not responsible for any injury or damage caused by use of this instrument for purposes other than those for which it is intended or by modifications of the instrument not performed by Bio-Rad Laboratories, Inc., or an authorized agent.

## Power Safety Information

### VOLTAGE SETTING INFORMATION

The universal hood of the Gel Doc XR+ or ChemiDoc XRS+ imaging system has a power supply that automatically chooses the correct voltage for your country or region.



## **FUSES**

The universal hood of the Gel Doc XR+ or ChemiDoc XRS+ imaging system has two user-serviceable fuses, F1 and F2, which are located on the bottom rear panel and are a part of the power entry module. Refer to the Maintenance chapter for fuse replacement instructions.







# Table of Contents

---

<b>Important Safety Information</b>	<b>i</b>
<b>Chapter 1. Introduction</b>	<b>1</b>
Gel Doc XR+ Imaging System	1
ChemiDoc XRS+ Imaging System	2
System Components	2
CCD Camera and Lenses	2
Universal Hood	3
Image Lab Software	3
Emission Filters	3
Optional Accessories	4
Printer	4
Conversion Screens	4
System Requirements	4
Computer Specifications	5
<b>Chapter 2. Gel Doc XR+ Imager</b>	<b>7</b>
Gel Doc XR+ Imager Workflow	7
Gel Doc XR+ Applications	8
Gel Doc XR+ Technical Specifications	9
<b>Chapter 3. ChemiDoc XRS+ Imager</b>	<b>11</b>
ChemiDoc XRS+ Imager Workflow	11
ChemiDoc XRS+ Applications	12
ChemiDoc XRS+ Technical Specifications	13
<b>Chapter 4. System Calibration</b>	<b>15</b>
<b>Chapter 5. Image Lab Software Overview</b>	<b>17</b>
Software Overview	18
Main Window	18
Toolbar	19
Results Data	19



Display Toolbox . . . . .	19
Start Page . . . . .	20
Analysis Tool Box . . . . .	20
Status Bar . . . . .	21
Menu Items . . . . .	21
<b>Chapter 6. Acquiring Images . . . . .</b>	<b>27</b>
Creating Custom Protocols . . . . .	27
Acquisition Settings . . . . .	28
Output Settings . . . . .	37
Choosing a Protocol . . . . .	37
Editing Protocols . . . . .	38
Position Gel . . . . .	39
Running a Protocol . . . . .	39
Application Tables . . . . .	40
<b>Chapter 7. Viewing Images . . . . .</b>	<b>43</b>
Results Overview . . . . .	43
Displaying Gel Images . . . . .	44
Display Gel Options . . . . .	44
Zoom Tools . . . . .	45
Fit in Window . . . . .	45
Image Transform . . . . .	46
Image Colors . . . . .	47
3-D Projection . . . . .	48
Image Info . . . . .	49
Displaying Data . . . . .	50
Lane and band definitions . . . . .	50
Volume definitions . . . . .	50
Analysis Table Options . . . . .	51
Lane Profile . . . . .	53
Standard Curve . . . . .	54
Report . . . . .	55
<b>Chapter 8. Analyzing Images . . . . .</b>	<b>57</b>
Analysis Tool Box Tools . . . . .	57
Auto Analysis . . . . .	57
Analysis Toolbox . . . . .	59
1. Image Tools . . . . .	60
2. Lane and Band Tools . . . . .	62



3. Molecular Weight Analysis Tools . . . . .	66
4. Quantity Tools . . . . .	67
5. Annotation Tools . . . . .	72
6. Volume Tools . . . . .	74
<b>Chapter 9. Generating Reports . . . . .</b>	<b>79</b>
Report . . . . .	79
Display Report General Tab . . . . .	80
Lane and Band Table Tab . . . . .	81
Volume Table Tab . . . . .	82
Print Report . . . . .	82
Print Report to .pdf File . . . . .	82
Adjust the Printer Settings . . . . .	82
<b>Chapter 10. Exporting Results . . . . .</b>	<b>83</b>
Exporting Gel Images . . . . .	83
Exporting Gel Images for Publication . . . . .	84
Exporting Gel Images for Analysis . . . . .	85
Exporting Gel Images to PulseNet . . . . .	85
Exporting Lane and Band Tables to Excel . . . . .	85
Exporting Volume Tables to File . . . . .	86
Snapshot Tool Export . . . . .	86
Analysis Table Export . . . . .	86
<b>Chapter 11. Maintenance . . . . .</b>	<b>87</b>
UV Transilluminator Lamp and Starter Replacement . . . . .	87
Epi-Illumination Lamp Replacement . . . . .	88
Fuse Replacement. . . . .	89
Calibrating Accessories. . . . .	93
Installing Optional Accessories. . . . .	93
UV/White Light Conversion Screen . . . . .	93
XcitaBlue™ Conversion Screen . . . . .	94
Gel Alignment Template Kit . . . . .	95
White Light Transilluminator . . . . .	97
Orange Fluorescence Reference Plate . . . . .	98
Ordering Information . . . . .	99
<b>Appendix A: Troubleshooting . . . . .</b>	<b>91</b>



**Appendix B: Accessories . . . . . 93**

Calibrating Accessories . . . . . 93

Installing Optional Accessories . . . . . 93

    UV/White Light Conversion Screen . . . . . 93

    XcitaBlue™ Conversion Screen . . . . . 94

    Gel Alignment Template Kit . . . . . 95

    White Light Transilluminator . . . . . 97

    Orange Fluorescence Reference Plate . . . . . 98

Ordering Information . . . . . 99

**Appendix C: Thermal Printer Settings . . . . . 103**

**Appendix D: Regression Calculation Methods . . . . . 107**

**Glossary . . . . . 109**

**Index . . . . . 111**



# 1 Introduction

---

Gel Doc™ XR+ and ChemiDoc™ XRS+ gel documentation systems offer high performance and are easy to use. They both contain a charge coupled device (CCD) camera to capture images in real time and allow you to accurately position your sample and generate optimized image data.

The Gel Doc XR+ and ChemiDoc XRS+ systems use the same light-tight enclosure (the universal hood), which contains built-in UV and white light illumination, but different CCD cameras. Both systems feature dynamic flat fielding technology for superior image uniformity and accurate quantitation.

Bio-Rad Image Lab™ software controls image capture and optimization for your selected applications, analyzes results, and produces reports based on your specified output, all in a single workflow.

## **Gel Doc XR+ Imaging System**

Gel Doc XR+ imaging system is a high-resolution gel documentation system that allows fast, easy quantitation of gels and blots. Position your sample inside the imager and follow the onscreen steps to run a protocol with only one click. You can customize your applications within an existing protocol or create a new protocol using many options presented by the Image Lab software.

Features include:

- Detection of dynamic ranges that cover > 3 orders of magnitude
- Dynamic flat fielding specific to fluorescent and colorimetric applications



- Smart, application-based protocol setup using Image Lab software, which assists you by presenting appropriate filter and illumination sources for each application
- Ability to upgrade to the ChemiDoc XRS+ system

## **ChemiDoc XRS+ Imaging System**

In addition to the gel documentation applications of Gel Doc XR+ presented above, the ChemiDoc XRS+ imaging system offers sensitive chemiluminescent detection. The system includes a supersensitive 16-bit CCD camera that is deeply cooled for faint sample detection and for accurate quantification of image data.

Features include:

- Imaging applications that require excellent sensitivity (chemiluminescent western blots)
- Smart, application-based protocol setup using Image Lab software, which assists you by presenting appropriate filter and illumination sources for each application
- Exceptional sensitivity and a dynamic range greater than 4 orders of magnitude
- Flexibility to image chemiluminescent, fluorescent, and colorimetric samples with dynamic flat fielding specific to each application

## **System Components**

### **CCD Camera and Lenses**

Either the Gel Doc XR+ or the ChemiDoc XRS+ camera is placed on top of a light-tight enclosure (the universal hood) for capturing images. The camera comes with a motorized zoom lens (MZL) that allows remote adjustment of the lens control functions (zoom, focus, and iris).

A patent-pending software algorithm controls the MZL, giving the user automatic image focus once an initial calibration is performed during system installation. See the Technical Specifications table, included in the chapter which discusses your camera, for complete specifications of each system.

A +1 diopter lens is factory installed to allow the entire sample stage to be visible. This lens should always remain on the MZL assembly.



## **Universal Hood**

The universal hood is designed to capture fluorescent and chemiluminescent images without using a photographic darkroom. The enclosure has built-in white light epi-illumination and UV transillumination. For easy sample loading, the UV transilluminator is located in the drawer of the universal hood and can be accessed from the front of the enclosure. When not imaging, the lights in the darkroom enclosure turn off automatically.

The universal hood has touchpad buttons to perform various functions; however, Image Lab software controls all of these functions remotely, removing any requirement for manual control of the lens and lights. Running a protocol overrides any touchpad input.

## **Image Lab Software**

Each system ships with a full version of Image Lab software. In addition to controlling the imager systems, image capture, and optimization, Image Lab software can be used to annotate and document images, analyze molecular weights (or base pairs, when imaging nucleic acid gels), and to determine accurate quantitation and purity of samples.

You can print all or a subset of your data in a report. Alternatively, you can export your data to other software, such as Microsoft Office programs, for further analysis or presentation options. See Chapters 5 to 10 for detailed instructions on the software.

## **Emission Filters**

The universal hood can hold two different emission filters for fluorescent applications. No filter is required to image chemiluminescent samples.

A standard filter (catalog # 170-8081) is used for colorimetric (white light) applications and is included in the installation kit.



## Optional Accessories

Bio-Rad Laboratories, Inc. offers a selection of optional filters and illumination sources. See Appendix B for ordering information and a complete listing of accessory filters, UV light sources, optional parts and replacement parts.

### Printer

For your convenience, Bio-Rad offers an optional USB printer for use with the Gel Doc XR+ and ChemiDoc XRS+ systems: the Mitsubishi P93DW Thermal Printer, (catalog #170-8089).

### Conversion Screens

#### WHITE LIGHT CONVERSION SCREEN

The white light conversion screen (catalog #170-8001) is a phosphor screen that produces white light transillumination when placed on top of the UV transilluminator.

#### WHITE LIGHT TRANSILLUMINATOR

An optional white light transilluminator (catalog #170-7950) may also be used with either system. The white light transilluminator plugs inside the universal hood and is controllable by the touchpad or the computer software.

#### XCITABLUE CONVERSION SCREEN

The optional XcitaBlue™ screen kit (catalog # 170-8182) is a UV to blue light conversion screen, which allows you to visualize DNA samples while protecting them from UV damage.

## System Requirements

Image Lab software runs under Windows XP Professional, Windows 7, and Mac OS X. Images scanned at high resolution can be quite large. The amount of memory required for using the program is mainly determined by the size of the images you will scan and analyze.

For this reason, we recommend that you archive images on a network file server or on removable storage media. Bio-Rad can also provide an appropriate computer to use with this system. Contact your local Bio-Rad representative for more details.



## Computer Specifications

Specifications	Minimum	Recommended
Operating system	Windows XP SP3	Windows XP SP3 Professional
	Windows 7, 32- and 64-bit	Windows 7 Professional, 64-bit
	Mac OS X 10.5 or 10.6	Mac OS X 10.6
Processor	Pentium 4 or equivalent (Windows) at 2.0 GHz	Core 2 Duo 2.0 GHz or higher
Hard disk space	20 GB	> 100 GB
Memory (RAM)	1,024 MB	> 1,024 MB
Ports for connecting instrument	1 free USB 2.0 port	1 free USB 2.0 port
Other software (optional)	Microsoft Excel 2000 or later (Windows)	Microsoft Excel 2000 or later (Windows)
	Office 2008 or iWork software (Mac)	Office 2008 or iWork software (Mac)







## 2 Gel Doc XR+ Imager

---

Refer to the Gel Doc XR+ Installation Guide (10016995) found in your Gel Doc XR+ Installation Kit (catalog # 170-8199) for instructions to install your Image Lab software and to assemble and calibrate your system.

To recalibrate your system because you have acquired new accessories, refer to Chapter 4.

### Gel Doc XR+ Imager Workflow

Following are the basic steps to acquiring, analyzing, and archiving an image using the Gel Doc™ XR+ system and Image Lab software:

1. Select an existing protocol or customize a new one.
2. Position the sample to be imaged.
3. Run your selected protocol.
4. View the displayed results.
5. Optimize the analysis.
6. Generate your report.
7. Save or export the results.



## **Gel Doc XR+ Applications**

The Gel Doc XR+ system is capable of running a wide variety of protocols. Those listed below are examples of the detection reagents that may be used to image your samples. Contact Bio-Rad technical support to determine whether your application can be imaged on this instrument.

See Chapter 6, Acquiring Images, for complete instructions on designing protocols.

### **NUCLEIC ACID GELS**

- Ethidium bromide
- SYBR® Green
- SYBR® Safe
- SYBR® Gold
- GelGreen™
- GelRed™
- Fast Blast™

### **PROTEIN GELS**

- Coomassie Blue
- Copper stain
- Zinc stain
- Flamingo
- Oriole
- Silver stain
- Coomassie Fluor Orange
- SYPRO Ruby
- Krypton

### **BLOTS**

- Colorimetric
- Qdots 525
- Qdots 565 Qdots 625
- CY2
- Alexa 488
- DyLight 488
- Fluorescein



# Gel Doc XR+ Technical Specifications

Applications	
Chemiluminescence	No
Fluorescence*	Yes
Colorimetry/densitometry	Yes
Gel documentation	Yes
Hardware Specifications	
Maximum sample size (cm)	28 x 36
Maximum image area (cm)	25 x 26
Excitation source	Trans-UV and Epi-White are standard (302 nm included, with 254 and 365 nm available as options). Optional Trans White, self-powered or conversion screen. Optional XcitaBlue™ UV/blue conversion screen.
Illumination control	3 modes Trans-UV, Trans White, Epi-White
Detector	CCD
CCD resolution (H x V)	1,392 x 1,040
Pixel size (H x V in microns)	4.65 x 4.65
Cooling system	Not applicable
Camera cooling temperature	Not available
Filter holder	3 positions (2 for emission filters, 1 for using no filter)
Emission filters	1 included (standard), 3 optional
Dynamic range	>3.0 orders of magnitude
Pixel density (gray levels)	4,096
Dynamic flat fielding	Application-specific, for all applications
Instrument size (cm)	36 (L) x 60 (W) x 96 (H)
Instrument weight (kg)	32
Operating Ranges	
Operating voltage	110/115/230 V AC nominal
Operating temperature	10–28°C (21°C recommended)
Operating humidity	<70% non-condensing
Automation Capabilities	
Workflow automated selection	Application driven, user selected or recalled by a protocol
Workflow automated execution	Controlled by a protocol via application specific setup for image area, illumination source, filter, analysis and reporting
Workflow reproducibility	100% repeatability via recallable protocols; from image capture to quantitative analysis and reports
Autofocus (patent pending)	Pre-calibrated focus for any zoom setting or sample height
Image flat fielding (patent pending)	Dynamic; pre-calibrated and optimized per application
Autoexposure	2 user-defined modes; (intense or faint bands)

\*Using the optional XcitaBlue kit (catalog #170-8182) is highly recommended performing SYBR® Safe DNA applications because the UV to blue conversion screen allows you to visualize DNA samples while protecting from UV damage.







## 3 ChemiDoc XRS+ Imager

---

Refer to the ChemiDoc XRS+ Installation Guide (10006999) found in your ChemiDoc Installation Kit (catalog # 170-8299) for instructions to install your Image Lab software and to assemble and calibrate your ChemiDoc XRS+ system.

To recalibrate your system because you have acquired new accessories, refer to Chapter 4.

To recalibrate your system because you have acquired new accessories, refer to System Calibration

### ChemiDoc XRS+ Imager Workflow

Following are the basic steps to acquiring, analyzing and archiving an image using the ChemiDoc™ XRS+ system and Image Lab software:

1. Select an existing protocol, or customize a new one.
2. Position the sample to be imaged.
3. Run your selected protocol.
4. View the displayed results.
5. Optimize the analysis.
6. Generate your report.
7. Save or export the results.



When running chemiluminescent blots, you can use a signal accumulation mode, which takes a series of exposures within a period of time. This enables you to choose an image with optimal signal. Refer to Chapter 6 Acquiring Images for instructions.

## ChemiDoc XRS+ Applications

The ChemiDoc XRS+ system is capable of running protocols to image all the detection reagents from the Gel Doc XR+ list, plus the others added to this list. Those listed are examples of the detection reagents that may be used to image your samples and not a complete list. Contact Bio-Rad technical support to determine whether your application can be imaged on this instrument.

See Chapter 6, Acquiring Images, for detailed instructions on designing protocols.

### NUCLEIC ACID GELS

- Ethidium bromide
- SYBR® Green
- SYBR® Safe
- SYBR® Gold
- GelGreen™
- GelRed™
- Fast Blast™ stain

### PROTEIN GELS

- Coomassie Blue
- Copper stain
- Zinc stain
- Flamingo
- Oriole
- SYPRO Ruby
- Silver stain
- Coomassie Fluor Orange
- Krypton

### BLOTS

- Chemiluminescent
- Colorimetric
- Qdots 525
- Qdots 565
- Fluorescein
- Qdots 625
- CY2
- Alexa 488
- DyLight 488



# ChemiDoc XRS+ Technical Specifications

Applications	
Chemiluminescence	Yes
Fluorescence*	Yes
Colorimetry/densitometry	Yes
Gel documentation	Yes
Hardware Specifications	
Maximum sample size (cm)	28 x 36
Maximum image area (cm)	25 x 26
Excitation source	Trans-UV and Epi-White are standard (302 nm included, with 254 and 365 nm available as options). Optional Trans White, self-powered or conversion screen. Optional XcitaBlue™ UV/blue conversion screen.
Illumination control	5 modes (Trans-UV, Epi-White, and No Illumination for chemiluminescence are standard) Trans White and XcitaBlue conversion screens are options
Detector	Supercooled CCD
CCD resolution (H x V)	1,392 x 1,040
Pixel size (H x V in microns)	6.45 x 6.45
Cooling system	Peltier cooled
Camera cooling temperature	–30°C controlled
Filter selector	3 positions (2 for filters, 1 without filter, for chemiluminescence)
Emission filters	1 included (standard), 3 optional
Dynamic range	>4.0 orders of magnitude
Pixel density (gray levels)	65,535
Dynamic flat fielding	Application-specific, for all applications
Instrument size (cm)	36 (L) x 60 (W) x 96 (H)
Instrument weight (kg)	32
Operating Ranges	
Operating voltage	110/115/230 V AC nominal
Operating temperature	10–28°C (21°C recommended)
Operating humidity	<70% non-condensing
Automation Capabilities	
Workflow automated selection	Application driven, user selected or recalled by a protocol
Workflow automated execution	Controlled by a protocol via application specific setup for image area, illumination source, filter, analysis and reporting
Workflow reproducibility	100% repeatability via recallable protocols; from image capture to quantitative analysis and reports
Autofocus (patent pending)	Pre-calibrated focus for any zoom setting or sample height
Image flat fielding (patent pending)	Dynamic; pre-calibrated and optimized per application
Autoexposure	2 user-defined modes; (intense or faint bands)
*Using the optional XcitaBlue kit (catalog # 170-8182) is highly recommended if performing SYBR® Safe DNA applications, because the UV to blue conversion screen allows you to visualize DNA samples while protecting against UV damage.	







## 4 System Calibration

---

When your system is installed, system calibration is performed with a calibration wizard. See the Installation Guide (10016995 for Gel Doc™ XR+ and 10016999 for ChemiDoc™ XRS+) in your Installation Kit for detailed instructions.

The instrument calibration wizard performs several procedures required to automate the system and prevent focus problems. Each of these calibrations affects your system as follows:

- **Focus Calibration**—This calibration allows automated focus settings at any zoom point, using a (patent pending) software algorithm. Therefore, your focus remains correct whether you view an entire sample or an area of interest.
- **Focus Calibration with Height Offset**—This calibration takes the tallest of the available conversion screens into account, and extrapolates values for the others, so that focus remains optimal, whichever screen is used.
- **Dark Reference Image** (Gel Doc XR+ only)—This calibration determines and corrects any background signal present in your imager.
- **Bias Reference Image** (ChemiDoc XRS+ only)—This calibration provides a bias value correction.
- **UV Flat Field Calibration**—This calibration generates the flat field correction profiles needed for the UV light source. Because of this calibration, your images have backgrounds of even intensity.
- **Lens Flat Field Calibration**—This calibration corrects for the intensity roll-off inherent in any lens.
- **White Conversion Screen Calibration**—This calibration generates a flat field correction profile needed for the white light conversion screen.



The Instrument Setup dialog box available from the Edit menu displays these calibration results for reference. The calibration wizard needs to be run only once, when your system is installed. The calibrations remain valid over the life of the instrument.

The calibration need not be changed unless you add equipment, such as a new light source.

When you add light sources or filters to a Gel Doc XR+ or ChemiDoc XRS+ system, recalibrate your system using the dialog box found at Edit > Instrument Setup.

**Instrument Setup**

**Hardware Information**

Model: ChemiDoc XRS+  
Serial Number: 721BR00019  
Camera Serial Number: A09H840004

**Illumination Options**

☒ White Light Conversion Screen ?  
☐ White Light Transilluminator ?  
☒ XcitaBlue Conversion Screen ?

**Filter Options**

Filter Position 1: Standard Filter ?  
Filter Position 2: No Filter ?

**Instrument Calibration**

**Focus Calibration:**  
Last Calibration: Thu Sep 17 12:37:50 2009 [Reset]  
Last Calibration (Offset): Fri Sep 18 13:50:00 2009

**Flat Field:**  
Last Calibration (Lens): Fri Sep 18 09:33:59 2009 [Reset]  
Last Calibration (White): Fri Sep 18 13:51:09 2009 [Reset]  
Last Calibration (Orange): Fri Sep 18 13:52:41 2009

**Bias Value Correction**  
Last Calibration: Thu Sep 17 12:38:04 2009 [Reset]

[OK]

1. If you have a new illumination source, select the appropriate box in the Illumination Options field.
2. If you are adding new filters, use the drop-down list to match what is installed in your instrument.
3. The software prompts you to reset the calibrations needed for the new illumination sources.
4. Click OK to exit the dialog box. Your settings remain until you make further changes.



## 5 Image Lab Software Overview

---

Image Lab™ image acquisition and analysis software works with the Gel Doc™ XR+ and ChemiDoc™ XRS+ imaging systems to create a reproducible, automated and time-saving workflow for running your gels.

In Image Lab software, a protocol is any combination of imaging, analysis, and report settings that has been saved to run as a single workflow. Researchers can run one protocol repeatedly or easily design a wide range of protocols.

With Image Lab software you can view analyzed data, edit the analysis, and produce customized reports that show precisely the settings applied, in order to ensure repeatable results.

Two types of files generated by Image Lab software:


- Protocol files, which describe the parameters for imaging and analyzing your gel images. Protocol files are saved with a .ptl extension
- Image files, which contain the imaged gel, annotations, and analysis performed on the gel. Image files are saved with a .scn extension

An imaged gel, run according to a protocol file, generates an image file.



## Software Overview

The following illustration shows the Image Lab software main window. The paragraphs below describe the main software elements.

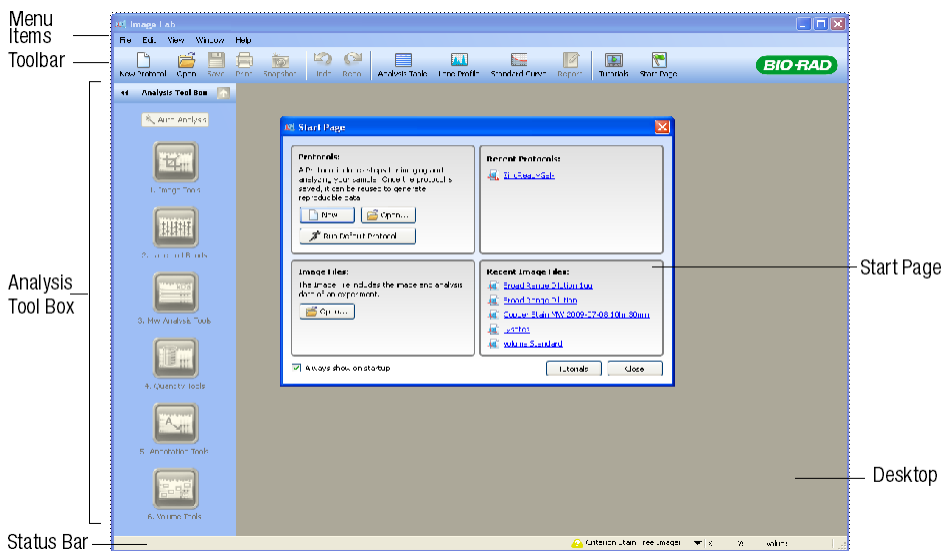

## Main Window

There is one main window used in Image Lab software. Image screens, protocol screens, and dialog boxes presenting choices all open on the desktop, the gray-colored area of the main window.

If you have many screens open on the desktop, you can make one visible by clicking on the dark blue bar at the top of the selected screen. A list of open protocols and/or image files is also found under the Window menu, and you can click your choice in that list to bring it to the top.

You can view all of the analysis for an image or protocol at once, or compare image results by arranging the screens on the desktop.




## Toolbar

Many tools available elsewhere in the software can be chosen directly by clicking on the toolbar icons. The Snapshot tool allows you to send a screen capture of your image to the clipboard, or save it as a file. You can view demonstrations of various functions by clicking on Tutorials. The unlimited Undo and Redo buttons allow you to correct missteps easily.

Clicking the green Default Protocols button at the right end of the toolbar displays the Default Protocol dialog box where you can set up, view, and edit default protocols.

Additional tools for viewing protocols and results data presented in various forms are labeled below. These tools are described in Viewing Images.

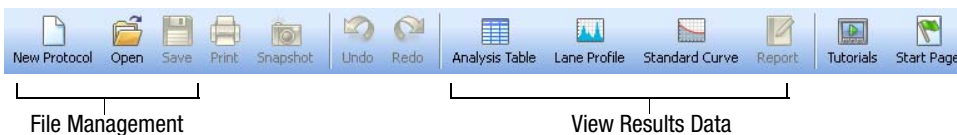


## Results Data

Results data associated with the gel images can be visualized several different ways; as an Analysis Table, a Lane Profile, a Standard Curve, or in a Report. These screens always display the analysis from the image you have selected.

Buttons to toggle these views on and off are located in the main toolbar, shown above. All of them can be viewed at one time. See “Displaying Data” on page 50 for details.

## Display Toolbox

The Display Toolbox near the top of every image allows you to display your images in the most useful ways. See “Displaying Gel Images” on page 44 for details.

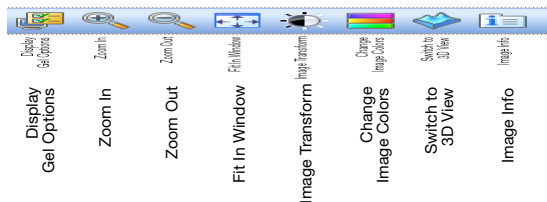



## Start Page

The Start Page guides users through creating, opening, and viewing protocols and images.



## Analysis Tool Box

The Auto Analysis button  quickly analyzes images.

The remaining tools customize the *analyzed* data.

**Note:** An image file must be selected (for example, if you are using a Windows computer, the title bar is dark blue) in order for any of these analysis tools to be available for use.

1. **Image Tools** allows the user to flip and rotate images, crop, and transform the image files.

2. **Lane and Band Tools** control the detection function, allowing the user to resize, adjust, and bend lanes, and detect, adjust, add or delete bands.

3. **MW (Molecular Weight)** (or base pairs, if using nucleic acid gels) analysis calculation allows the user to choose standard samples, assign standard lanes, and choose a regression method.


4. **Quantity Tools** allow the user to automatically quantitate bands, using either relative or absolute values.



5. **Annotation Tools** are useful for drawing attention to any area of a gel that is of interest to the researcher.
6. **Volume Tools** allow you to manually quantify an object inside a boundary that you define.

These tools are described in “Analysis Tool Box Tools” on page 57.

## Status Bar

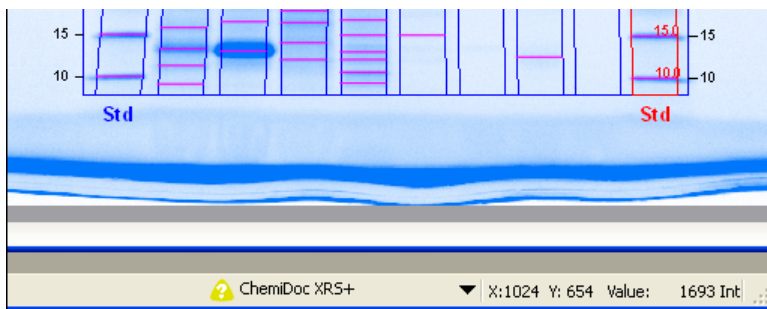

The status bar in the lower right-hand corner of the main window shows the imager that is being used, and the X and Y values for the cursor position on an image file. The Int (intensity) values show a numerical value for the image intensity at the mouse position. The maximum data range using a Gel Doc XR+ system is 0–4,095 (ChemiDoc XRS+ maximum data range is 0–65,535) but the ranges will vary according to the values contained within your image.

## Menu Items

The following section describes all of the choices under the Menu Items: File, Edit, View, Window, and Help. Many of these functions are also available in other locations, such as the Toolbar or the Start Page.

### FILE

File > **New Protocol** allows the user to create a new protocol through one window which contains all of the necessary steps and choices. See the Protocols chapter for detailed instructions. The Protocol can also be altered and stored for reuse.

File > **Open** browses the user's file system to retrieve a previously-saved protocol file or image file.

File > **Recent Images** gives the user a choice of opening a recent image file.

File > **Recent Protocols** gives the user a choice of opening a recent protocol.



File > **Save** allows the user to save a protocol or image file in their file system, once it is named.

File > **Save As** allows the user to name and store a protocol or image in their file system. Protocols are stored with a .ptl file extension. Image files are stored with a .scn extension.

File > **Close** gives the user a choice of closing the active desktop screen.

File > **Close All** closes all of the desktop screens.

File > **Export** enables the user to export gel images or analysis tables with the following options:

- **Export for Publication**—Exports a displayed image to a file. You can select from .bmp, .png, .jpg, and TIFF formats. The gel displays with any lanes, bands, and annotations that appear on the screen. See Exporting Results for more information.
- **Export for Analysis**—Creates a TIFF formatted file that retains all gel image data. Analysis data are not included. Use this option to analyze the image in other software such as Quantity One<sup>®</sup>, FPQuest<sup>™</sup>, or InfoQuest<sup>™</sup>FP.
- **Export for PulseNet**—Reduces the image to an 8-bit TIFF image file. Resolution is limited and file size is restricted to 300 dpi.
- **Lane and Band Table to Excel**—If you have Excel installed on your computer, it will launch with your Lane and Band Table displayed as a spreadsheet.
- **Lane and Band Table to File**—This choice exports as a CSV (comma-separated values) file, so your lane and band table can be opened in a database application.
- **Volume Table to Excel**—If you have Excel (or Numbers, on a Mac) installed on your computer, it will launch with your Volume Table displayed as a spreadsheet.
- **Volume Table to File**—This choice exports as a CSV (comma-separated values) file, so your volume table can be opened in a database application.

See Chapter 10 for more information about exporting files.



File > **Image Info** shows information about the individual gel and blot images, such as acquisition date, the data range, and image capture details, such as the exposure time and illumination source used. There are three tabs (Image Details, Analysis Settings, and Notes) which reveal specific information.

The screenshot shows the 'Image Info - Blot' dialog box with the 'Image Details' tab selected. The dialog is divided into two main sections: 'Acquisition Information' and 'Image Information'. Each section contains several text input fields with pre-filled values.

Acquisition Information	
Imager	ChemIDoc XRS+
Exposure Time (sec)	0.317 (Auto - Intense Bands)
Flat Field	Applied (Lens)
Serial Number	721BR00019
Software Version	2.0
Application	Colorimetric
Excitation Source	White epi illumination
Emission Filter	Standard Filter

Image Information	
Acquisition Date	9/18/2009 11:35:49 AM
User Name	Researcher
Image Area (mm)	X: 95.0 Y: 71.0
Image Pixels	X: 1392 Y: 1040
Pixel Size (um)	X: 68.2 Y: 68.2
Data Range (Int)	13665 - 64753

At the bottom of the dialog are 'OK' and 'Cancel' buttons.

File > **Page Setup** contains print controls, such as orientation (landscape or portrait), margins, printer used, and paper size.

File > **Print** displays a print preview of the gel and the header information, which includes the filename of the image, the user's name, and date/time it was printed. The usual Windows Print screen is available as well, which lets you select a printer, and the number of copies.

File > **Exit** closes Image Lab software (after asking you to save changes on your protocols or images).



## EDIT

Edit > **Undo** undoes the last user action.

Edit > **Redo** restores the last user action, after an Undo.

Edit > **Snapshot** allows the user to take a screen capture of the Lane Profile Window, the Standard Curve Window, or the default choice, Current Image View. This capture can include the name of the image, and it can be placed on the clipboard, or saved in a file.

Edit > **Default Imager** allow users who own two or more imagers to switch between them.

Edit > **Instrument Setup** allows users to review the instrument serial number and how their imaging system is calibrated. If accessories are added, use the Instrument Setup dialog box to change the system calibration to use them.

Edit > **Report Settings** allows you to configure your reports. This dialog box contains three tabs. All of the checkboxes are selected by default; clear boxes for information you do not want to include in reports. Your selections apply to all reports, until you change them again.

- The General tab has options for excluding or reporting information about your gel image.
- The Lane and Band Table tab allows the researcher to choose whether to include all, or just some lanes, with appropriate identifiers. Lane profiles can also be included.
- The Volume Table tab allow the researcher to choose appropriate identifiers for the Volume Table, and to exclude the table from his reports, if desired.

Edit > **Preferences** contains two tabs.

- The Protocol tab contains some presets for naming your image files. You can choose to include a designated Prefix, User Name, Date and/or Time in the name of your image files.
- The Colors tab allows you to choose colors for the graphic elements in your gels, such as Lane Frame, Lane, Band, Band Attribute and MW Legend. This functionality ensures that these elements are visible, whatever colors your gels are.



## VIEW

View > **Image Overview** opens a small window which shows your entire gel image, and a red rectangle outlining the area visible in the larger main window. This is useful when you are zoomed into a small section of your image.

View > **Image Transform** opens a window showing a histogram, where you can adjust the light and dark values of your gel image. This does not change your data, only the way it displays on your monitor.

View > **Operations History** opens a window showing sequential actions performed both by the user and the software.

## WINDOW

The Window controls allow you to show and hide multiple open image files on your desktop. A list of all currently open images and protocols display in this menu.

Window > **Tile** aligns all open image files so that they are all visible at once.

Window > **Tile Horizontal** places all open image files from top to bottom.

Window > **Tile Vertical** places all open image files from left to right.

Window > **Cascade** stacks all open image files and protocols with overlapping title bars, so that each one can be easily chosen for view.

Window > **Imitate Zoom** changes the zoom setting of all open images to the same zoom as the current image file.

Window > **Imitate Transform** changes the brightness and contrast of all open images to the same transform settings as the current image file.

Window > **Next** cycles through all open image files from oldest to newest.

Window > **Previous** cycles through all open image files from newest to oldest.

## HELP

Help > **Image Lab Help** displays online help.

Help > **User Guide** displays the user manual, available in .pdf form.

Help > **About** displays Image Lab software version and release date.







## 6 Acquiring Images

---

Image Lab™ software can be used to run specific applications with repeatable workflows, and to design custom protocols with a wide variety of settings that can be retrieved, revised, and reused.

In Image Lab software, a protocol is any combination of settings for imaging, analyzing and reporting that has been saved to run as a single workflow.

### Creating Custom Protocols

You start a new protocol by choosing the New Protocol box on the Start Page, or by choosing New Protocol from the toolbar or menu bar. Creating custom protocols involves choosing parameter settings from each screen.

#### PROTOCOL SUMMARY

The Protocol Summary screen can be used to quickly review all protocol settings.

#### PROTOCOL CHOICES

Protocols involve choices under three major processes: Choosing Acquisition Settings, Analyzing the Image, and Choosing Report Settings.

There are steps under each process where parameters are set. The choices available under each step appear as each Protocol Step is chosen.


If you want to make choices under a protocol step, click its description, for example, “3. Analyze Molecular Weight”.

To enable or disable any step, click its checkbox.




## Acquisition Settings

### STEP 1. GEL IMAGING

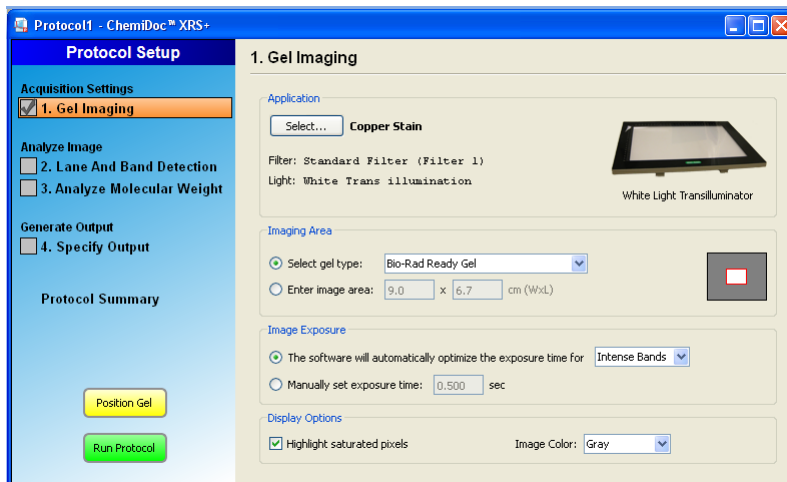

### APPLICATION OPTIONS

1. **Choose an application** from a menu that displays when you click the Select button.

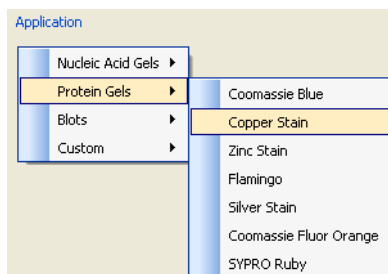


The applications are listed in submenus, under four application types. When you choose your application, any required filter or illumination source displays on the Protocol Setup screen.


Use the Custom application either when you want to run an existing application with a new name, or when you have an application unlike any existing applications. Selecting Custom from the applications menu displays the Manage Custom Applications dialog box. If you have stored Custom Applications, they display here.



To create a new custom application, select New from the Manage Custom Applications dialog box.

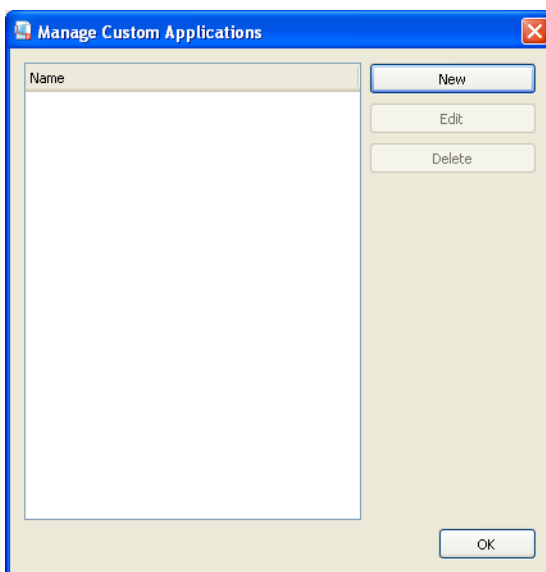


The Create Custom Applications dialog box appears.

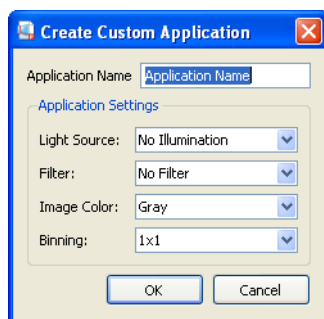

2. Choose a unique application name.
3. Select a light source, filter, and image color from the dropdown menus.
4. ChemiDoc XRS+ only) Select a binning setting. . Choosing a higher binning setting combines pixels to increase the amount of signal without increasing noise. While a higher setting provides optimal sensitivity for low-light applications such as chemiluminescence, it also reduces image resolution.




5. Click OK to create your custom application.
6. **Choose the Imaging Area** from a list of Bio-Rad gels, or enter image area dimensions. The red line represents the imaging area for the selected gel, and the gray rectangle represents the imager sample stage.

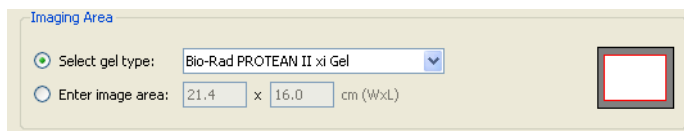


The 'Imaging Area' dialog box contains two radio buttons. The first, 'Select gel type:', is selected and points to a dropdown menu showing 'Bio-Rad PROTEAN II xi Gel'. The second, 'Enter image area:', is unselected and points to two input fields containing '21.4' and '16.0' followed by 'cm (WxL)'. To the right is a diagram of a rectangular sample stage with a red rectangle inside representing the imaging area.

7. For a list of applications with all required detection reagents, light sources, and any conversion screens or filters noted, see page 40. **Choose Image Exposure Time**—If you select the top radio button, Image Lab software acquires an image estimating an optimal exposure time. This option ensures the best use of the dynamic range.
  - If you choose Intense Bands from the dropdown menu, exposure is optimized for all bands.
  - Choosing Faint Bands from the dropdown menu results in a longer exposure where faint bands are more visible, but more prominent bands may be overexposed.
  - Select the manual exposure radio button to override automated imaging. That may cause saturation (overexposure) of more prominent bands. Exposure time can range from 0.001 to 99 seconds using the Gel Doc XR+, and from 0.001 to 7200 seconds, using the ChemiDoc XRS+.

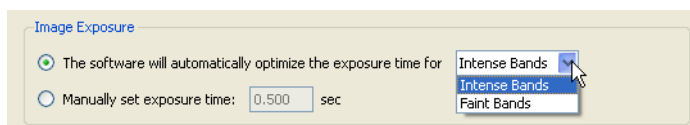

The 'Image Exposure' dialog box has two radio buttons. The first, 'The software will automatically optimize the exposure time for', is selected. It points to a dropdown menu currently showing 'Intense Bands'. The second radio button, 'Manually set exposure time:', is unselected and points to an input field showing '0.500' followed by 'sec'.

After imaging a gel with automatic exposure optimization, the used exposure time displays in the protocol (see below), so that you can manually adjust the exposure time if needed.

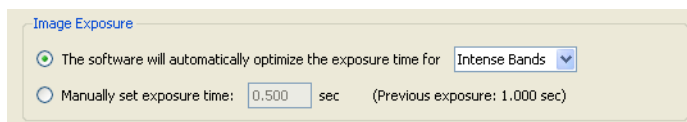

The 'Image Exposure' dialog box is shown after an exposure. The first radio button remains selected and points to the 'Intense Bands' dropdown. The second radio button, 'Manually set exposure time:', is now selected and points to an input field showing '0.500' followed by 'sec'. To the right of this field, text indicates '(Previous exposure: 1.000 sec)'.

You can also view the exposure time of the image later, in the Image Info window (refer to “Image Info” on page 49).




**Signal Accumulation Mode**—If you are running a chemiluminescent application on the ChemiDoc XRS+ system, you can also use signal accumulation mode (SAM). SAM is used to simplify capture of a good image of a chemiluminescent sample. This sample type often requires long integration times to obtain an image that represents the best range of signal.

Rather than manually acquiring a series of independent images with different imaging times, SAM presents a series of cumulative images with progressively greater signal in each image. SAM requires the user to estimate the shortest and longest times expected to generate an image with desired signal intensity. The user then decides how many total images they wish to acquire in this window of time.

For example, if the minimum time expected to image the sample is one minute and the maximum is 5 minutes, these values are entered (in seconds) in the setup window. By entering 5 in the Total number of images field, 3 images will be acquired between the first and last images.

The bar in the dialog box below shows that images will be acquired at one minute intervals, starting at one minute and ending at five minutes. The second one minute image is added to the first one minute image, resulting in a two minute total integration time image. The third one minute image is added to the previous image, and so on, until the last image is presented.

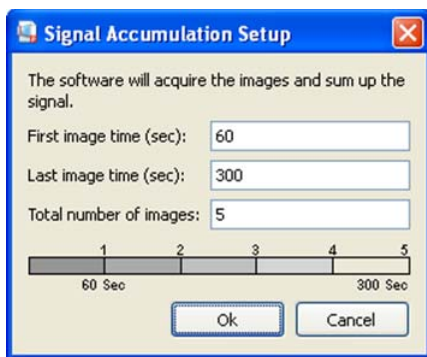

The software will acquire the images and sum up the signal.

First image time (sec):

Last image time (sec):


Total number of images:

1 2 3 4 5  
60 Sec 300 Sec

Ok Cancel



The most recent image is displayed by default in the main window.

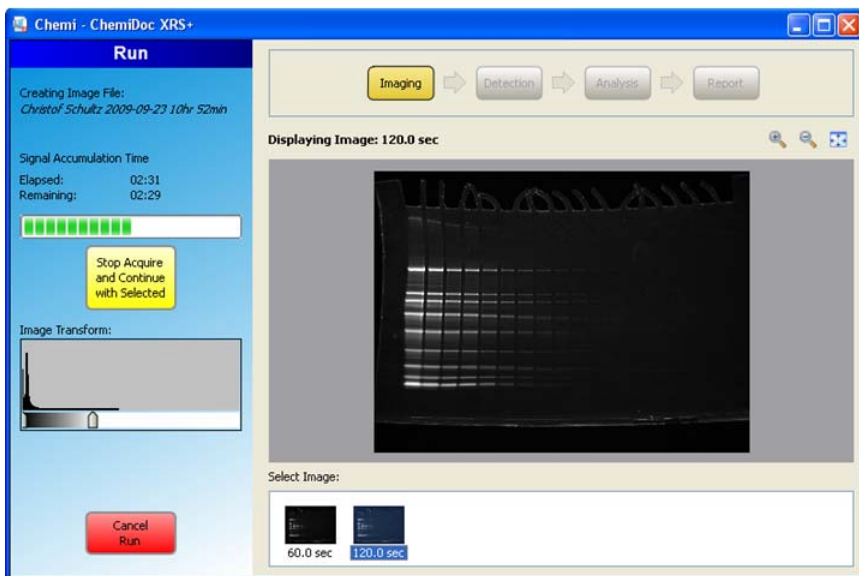

You can cancel the run by clicking on the red “Cancel Run” button, which ends the protocol. An image can be saved and analyzed at any time by right-clicking on the thumbnail and selecting Save from the pop up menu. Then click the yellow button labeled “Stop Acquire and Continue with Selected”. This choice stops the acquisition and continues the protocol using the selected image. All other images acquired will be discarded.

Although SAM is useful to determine the optimum imaging time for a chemiluminescent sample, it does not result in data as accurate as a single image. Signal that is near the intensity of background noise becomes increasingly masked as the number of cumulative images grows. Therefore, if you need to identify extremely faint signals in your image, reacquire it as a single image, using the time the SAM tool found to be appropriate.

## 8. Choose Display Options


**Highlight saturated pixels**—Select the Highlight saturated pixels checkbox to see any saturated pixels in red, in order to show how much of the gel image is saturated. This option can be changed later by selecting View > Image Transform.

**Image Color**—Select color choice to display the sample image. Viewing the image with a different color scheme can make it easier to see all of the elements. See “Image Colors” on page 47 to see the color choices dialog box.



## Analyze Image

### STEP 2. DETECT LANES AND BANDS

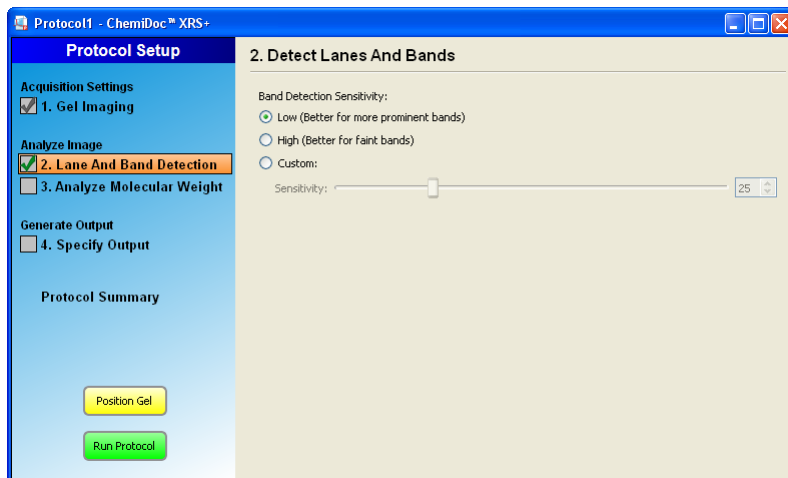

To perform any analysis of the gel or blot, Image Lab software needs to detect lanes and bands on the image. Lanes are detected automatically, and then the background is auto-subtracted. Refer to Lane and Band Tools on page 62 for full details. You have the following choices to customize band detection:

#### LOW BAND DETECTION SENSITIVITY

This setting will set detection at a low level, for images with more prominent bands. Faint bands will not be detected with this setting.

#### HIGH BAND DETECTION SENSITIVITY

This setting will set detection at a higher level, for images that are more faint. Extraneous bands can be removed later, using the Band Tools in the Analysis Tool Box. See “2. Lane and Band Tools” on page 62.

#### CUSTOM

Select a numerical value between 1-100, in order to choose the best detection sensitivity for your sample.


**Note:** When Low or High Band Detection Sensitivity is used, these are the numerical values that are set:

Low sensitivity = 25



High sensitivity = 75

### STEP 3. ANALYZE MOLECULAR WEIGHT

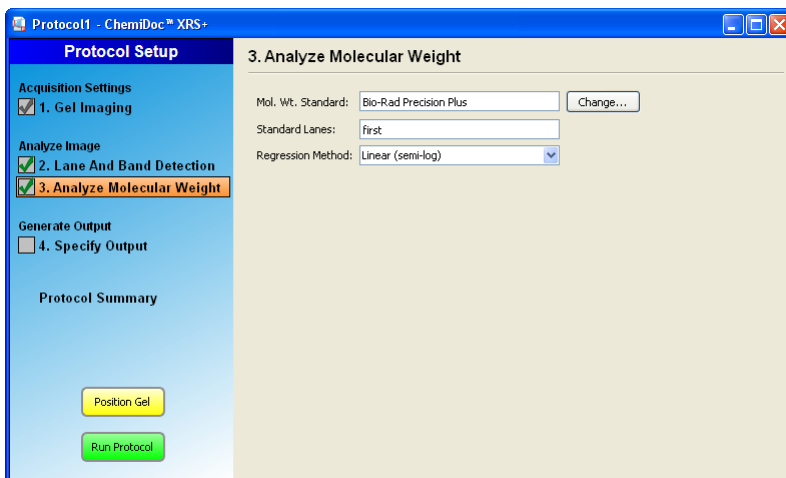

To estimate the size of the molecules in the bands of your gel, use this step to enter the standards you are using, and also to specify the lane(s) in which the standards are placed. For nucleic acid samples, use this step to determine the size of the bands in base pairs.

You can analyze molecular weight automatically by clicking Protocol Step 3. “Analyze Molecular Weight”. If this box is selected, the software calculates the molecular weight (or base pairs) for each band, based on the specified standard.

### STANDARDS

Determining molecular weight depends upon the selection of the proper protein standards. Many protein standards are available from Bio-Rad including the following:


- Precision Plus Protein™ Standards
- Prestained SDS-PAGE Standards Broad Range
- Prestained SDS-PAGE Standards Low Range
- Prestained SDS-PAGE Standards High Range



Many different DNA standards are also available. See the Edit Button illustration below for a partial list of DNA standards. See the Ordering Information on page 99 for all of the standards available from Bio-Rad and their part numbers.

### EDIT BUTTON

The Edit button opens the Manage Standards dialog box, which allows you to add other third-party standard samples, and to edit your list.

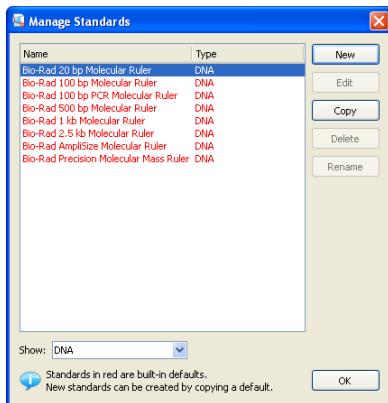

### STANDARD LANES

You can choose which lane contains your standards by typing lane numbers or the words First and Last in the Standard Lanes field. The format is xx, xx, xx, ... where xx is the lane number. For example, if you run an 18-well gel and want your standards in lanes 1, 10 and 18, enter First, 10, and Last.

**Note:** Lane detection works best when standards are placed in the first and last lanes.



## REGRESSION METHODS

A regression method is used to calculate the molecular weight/size of the unknown bands. The software will use the relative front and molecular weight values of the standard bands to calculate the standard curve. This standard curve is then used to calculate the values of the unknown bands. The shape of the standard curve is based on the selected regression method. You have a choice of four regression methods.

The methods and the minimum number of standard bands required for each type are shown in the following table.

Regression Method	Minimum number of standard bands
Linear (semi-log)	2
Point-to-Point (semi-log)	2
Logistic	5
Cubic Spline	5

If you do not have enough data points for the selected method, the molecular weight of the unknown bands will not be calculated.

Which method to select depends on the gel type:

- **Gradient gels:** The Linear (semi-log) regression method works well for these gels, since the mobility of the bands is linear to the log of their molecular weight. As an alternative, the Point-to-Point (semi-log) method can be used, if the  $R^2$  value is not sufficient.
- **Fixed percentage gels:** These gels have a non-linear relationship between the mobility and the molecular weight. Good choices for regression methods are Logistic or Cubic Spline.

You can check how well each regression method fits the data in the standard curve window (for more information, see “Standard Curve” on page 54). The Linear (semilog) regression method provides a measurement which describes how well the standard curve fits the data,  $R^2$  value. The closer the  $R^2$  value is to 1.0, the better the data fits the standard curve.

The molecular weight/size of each band is displayed in the Analysis Table in the Mol. Wt./Base Pair column. Refer to page 67.

For information about the calculations behind the regression methods, see “Appendix D: Regression Calculation Methods” on page 107.



## Output Settings

Image Lab software prints to the default printer, unless you select otherwise.


You have three choices for specifying output:

- You can automatically print the image; this is useful for gel documentation, especially if a thermal printer is connected. See Appendix C: Mitsubishi P93/P95 Thermal Printer Settings, for instructions on setting up the Mitsubishi P93/P95 thermal printer.
- You can automatically print the report, including all accompanying analysis information.

**Note:** You cannot print a report on a thermal printer.

- You can display the image on your computer.

### STEP 4. SPECIFY OUTPUT

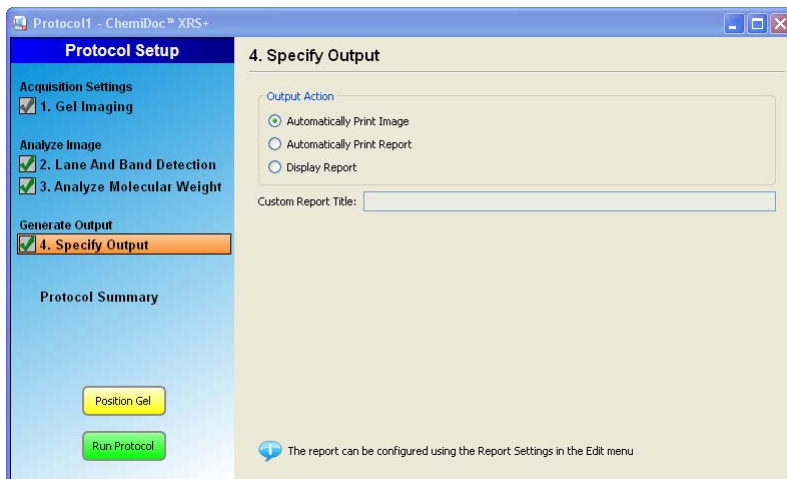

Here you can choose whether to automatically display or print your report after your protocol ends.

The report can be customized with the options in Edit > Report Settings.

For information about reporting options, see “Report” on page 79.

## Choosing a Protocol

Use the same protocol repeatedly by clicking the Open button on the Start Page, or by clicking File > Open from the menu bar.




You can also choose recently used protocols or image files from the lists on the Start Page.



## Editing Protocols

Existing protocols can be changed and/or renamed using the many tools in Image Lab software.


Open any protocol you have saved in the past by choosing File > Open. You will be presented with the same set of menus and choices described in the previous section, “Creating Custom Protocols” on page 27.

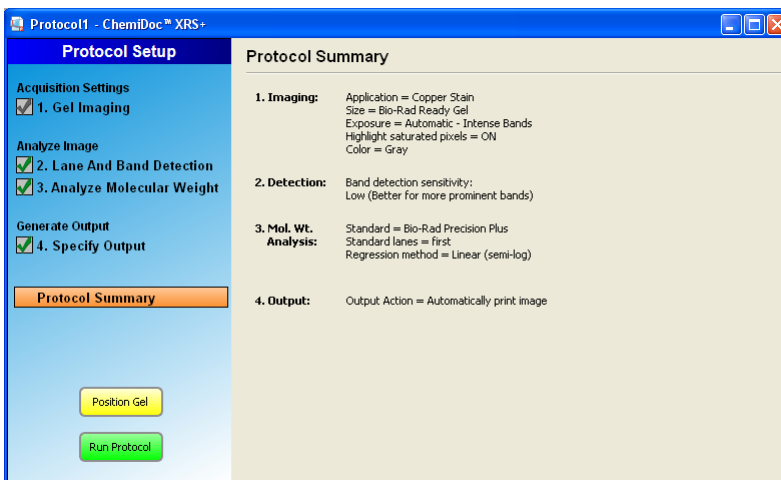

Make and save your changes without renaming the protocol to replace an older one. Use “Save As” with a different name to make a second protocol.



## Position Gel

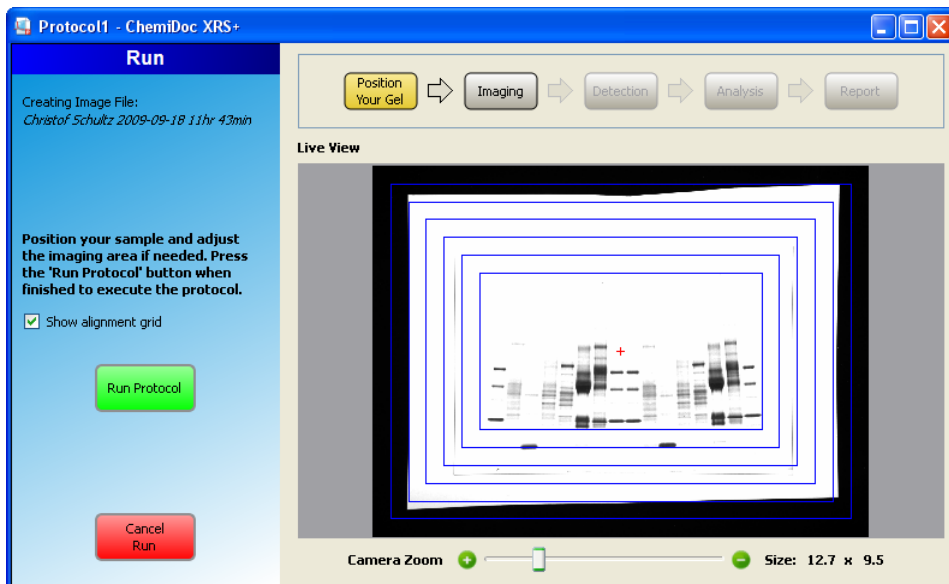

To position a gel, perform these steps:

1. Click the yellow Position gel button from the Protocol Setup field shown on page 38.
2. Place your gel on the imager stage and view your gel within Image Lab software.
3. Use the slider below the image to zoom your image into place. You may also move your gel manually until it is centered properly on the stage.

**Note:** The Bio-Rad gel alignment template kit supports four sizes of standard agarose gels to be centered quickly and easily. Refer to page 95 in the Accessories chapter for more information.

## Running a Protocol

To execute a protocol, click the green Run Protocol button shown above.



## Application Tables

The following tables contain all of the light sources, conversion screens, or filters you need for each application.

**Table 1. Nucleic acid gels**

Detection Reagent	Preferred Light Source	Alternate Light Source	Preferred Filter	Alternate Filter	Preferred Flat Field	Alternate Flat Field
Ethidium bromide	UV	None	Standard	630/30	Orange	Lens
SYBR® Green	XcitaBlue™	UV	560/50	Standard	Lens	None
SYBR® Safe	XcitaBlue	UV	560/50	Standard	Lens	None
SYBR® Gold	XcitaBlue	UV	560/50	Standard	Lens	None
GelGreen	XcitaBlue	UV	560/50	Standard	Orange	Lens
GelRed	UV	XcitaBlue	630/30	Standard	Lens	None
Fast Blast™ stain	White Trans	None	Standard	None	White Conv	Lens

**Table 2. Protein gels**

Detection Reagent	Preferred Light Source	Alternative Light Source	Preferred Filter	Alternative Filter	Preferred Flat Field	Alternate Flat Field
Coomassie Blue	White Trans	None	Standard	None	White Conv	Lens
Copper stain	White Trans	None	Standard	None	White Conv	Lens
Zinc stain	White Trans	None	Standard	None	White Conv	Lens
Flamingo	UV	None	Standard	Standard	Orange	Lens
Oriole	UV	None	Standard	None	Orange	Lens
Silver stain	White Trans	None	Standard	None	White Conv	Lens
Coomassie Fluor Orange	XcitaBlue	UV	560/50	None	Orange	Lens
SYPRO Ruby	UV	None	Standard	None	Orange	Lens
Krypton	XcitaBlue	UV	560/50	Standard	Orange	Lens



**Table 3. Blots**

<b>Detection Reagent</b>	<b>Preferred Light Source</b>	<b>Alternative Light Source</b>	<b>Preferred Filter</b>	<b>Alternative Filter</b>	<b>Preferred Flat Field</b>	<b>Alternate Flat Field</b>
Chemiluminescent* (only on ChemiDoc XRS+)	None	None	None	None	Lens	None
Colorimetric	White Epi-	None	Standard	None	Lens	None
Qdots 525	UV	None	520/30	Standard	Lens	None
Qdots 565	UV	None	560/50	Standard	Lens	None
Qdots 625	UV	None	630/30	Standard	Lens	None
CY2	UV	None	520/30	Standard	Lens	None
Alexa 488	UV	UV	520/30	Standard	Lens	None
DyLight 488	UV	UV	520/30	Standard	Lens	None
Fluorescein	UV	None	520/30	Standard	Lens	None








## 7 Viewing Images

Once your gel has been imaged, the image appears on the desktop. There are many controls to optimize viewing and to analyze your image.

### Results Overview

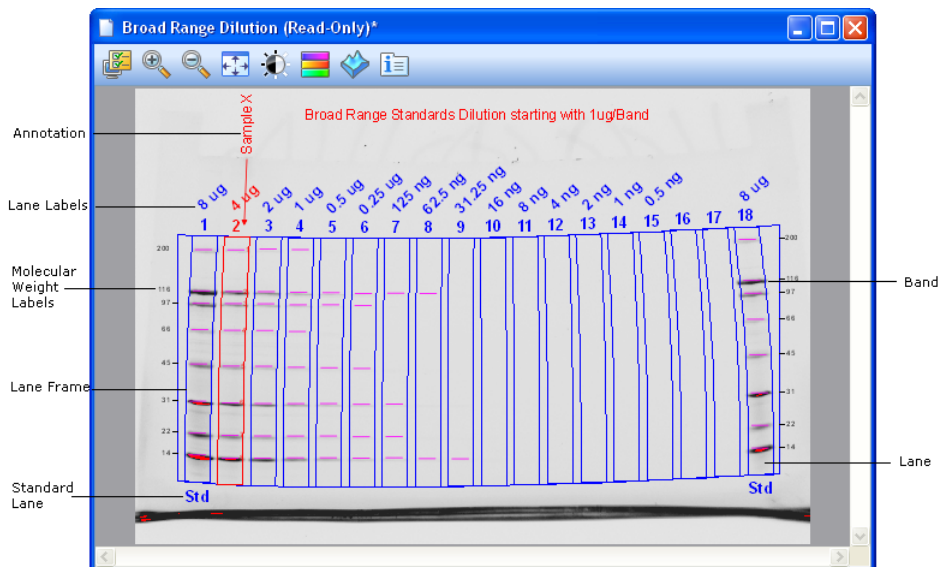

The above image shows a gel image with band and lane detection, as well as annotations. The labels show overlays that you can show or hide. Refer to “Display Gel Options” on page 44 for more information.




There are many ways to view the data associated with your results. You can view your data as an analysis table, a lane profile, a standard curve and as a report. See “Displaying Data” on page 50.

## Displaying Gel Images

Notice the display toolbar icons above the gel image on page 39. Each of these tools are described in the following sections.

### Display Gel Options

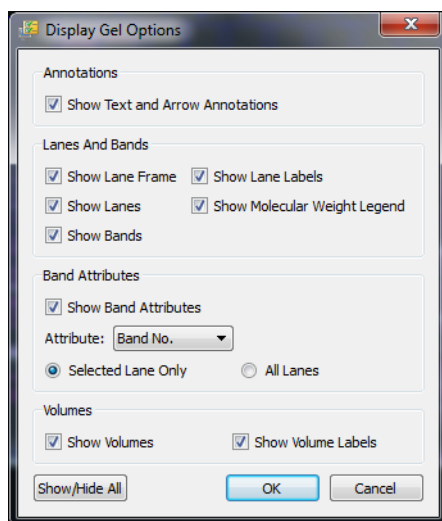

#### ANNOTATIONS

You have a choice whether or not to show text and arrow annotations that have been drawn on the image.

#### LANES AND BANDS

You can turn any image overlays, such as lane frames, lanes, bands, lane labels, and molecular weight legends on or off.



## BAND ATTRIBUTES

You can show the following attributes for selected lanes or for all lanes.

- Band number
- Band label
- Molecular weight
- Relative front
- Volume
- Abs. Quant.
- Rel. Quant.
- Band %
- Lane %

## VOLUMES

If you have drawn volume boundaries on your gel, you can display the volume boundaries and their volume labels.

## Zoom Tools



The zoom tools resize your gel image. Click on the magnifying glass with the plus to make your image larger; the minus to make your image smaller.

You can also zoom with the right mouse button. Click and drag the right mouse button to select an area you want magnified. Click the right mouse button again to return to the original view.

You can also resize your image with the scroll wheel on your mouse, if your mouse has one.


## Fit in Window



If you are currently zoomed in on an area of your image, clicking this button will bring your entire image into view.



## Image Transform

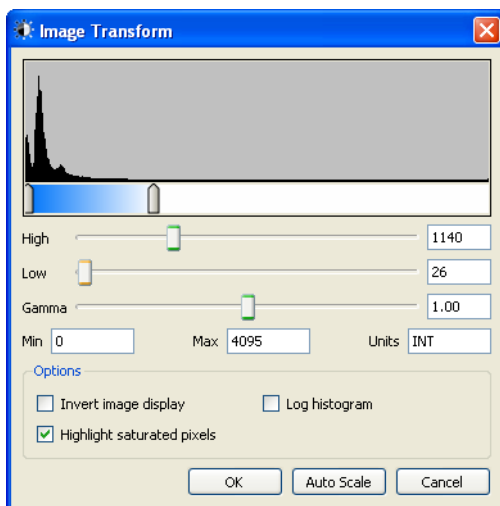

Use the Image Transform dialog box to adjust image brightness and contrast, optimizing the image display so faint detail can be seen. The minimum to maximum range will vary, according to the light and dark values present in your image.

**Note:** These adjustments do not change the data, only the way the data are displayed. The human eye cannot see as much range as is contained in the image.

The frequency distribution histogram shows the total data range in the image and the amount of data at each point in the range.

The Auto Scale button will determine an optimal setting for your particular image automatically. The lightest part of the image is set to the minimum intensity, and the darkest is set to the maximum.

- The High slider determines which intensity value will be shown at the maximum grayscale (or other color) in your gel image
- The Low slider determines which intensity value will be shown at the minimum grayscale (or other color) in your gel image
- The Gamma slider changes the grayscale curve. A value of 1 is linear. A value less than 1 redistributes a greater proportion of the grayscale to the first half of the intensity values. A value greater than 1 redistributes a greater proportion of the grayscale to the second half of the intensity values.




You can also type numerical values in the text boxes next to the sliders. Clicking anywhere on the slider bars will move the sliders incrementally.

## Options:

- **Invert image display**—inverts dark bands on a light background to light bands on a dark background, and vice versa
- **Highlight saturated pixels**—when this checkbox is selected, areas of the image with saturated (above a measurable range) signal intensity are highlighted in red
- **Linear or Logarithmic histogram**—This adjustment changes the Y axis on the histogram to display how many pixels are at each intensity value

## Image Colors

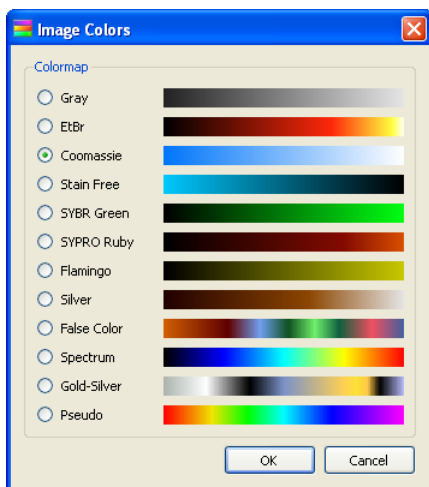


This dialog box allows you to choose a colormap for your image results file. Viewing the image with a different color scheme can make it easier to see all of the elements in the image, but does not change your data.



You have many color choices. The first eight imitate the colors of stained gels. The remaining choices supply enough color variation to highlight small differences in the image data.

- Gray
- EtBr (Ethidium bromide)
- Coomassie
- Stain Free
- SYBR Green
- SYPRO Ruby
- Flamingo
- Silver
- False Color
- Spectrum
- Gold Silver
- Pseudo

### 3-D Projection

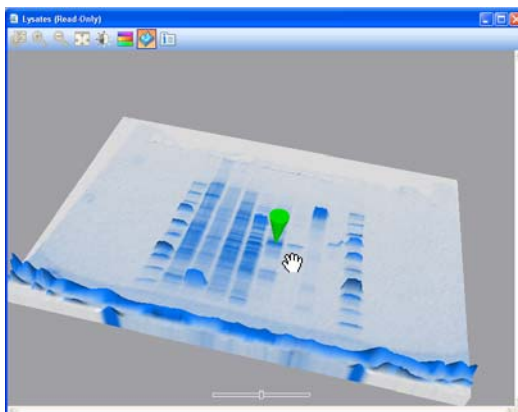

Clicking the 3-D icon transforms your gel image to a solid 3-dimensional model spinning in space with x, y and z dimensions. Accentuate or diminish the relative heights of data points by pulling the slider at the bottom of the window to the right or left.

To open the viewer, select the 3-D icon from the toolbar above your image.

Use these steps to view the intensity of various bands:

1. Click and drag the model to rotate into your preferred view.
2. Bring the window into focus by clicking on the image.
3. Press C to display a green cone, which can be dragged around with your cursor to evaluate the intensity of various bands.



## Image Info

The Image Info dialog box provides information about an active image. Click the icon on the image toolbar to display three tabs of information about your image.

### IMAGE DETAILS

Acquisition and image information is displayed in this dialog box.

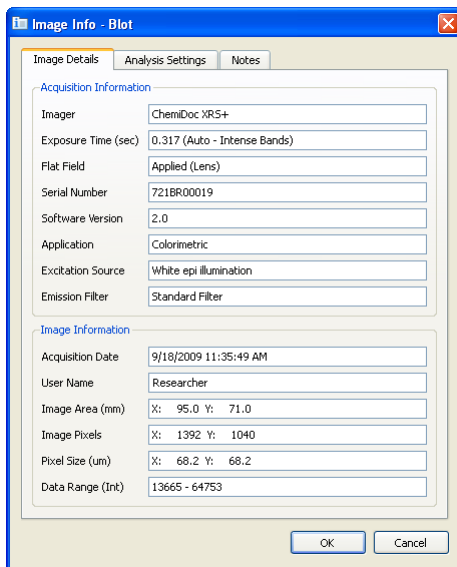


Image Info - Blot	
<div> <div>Image Details</div> <div>Analysis Settings</div> <div>Notes</div> </div>	
Acquisition Information	
Imager	ChemIDoc XRS+
Exposure Time (sec)	0.317 (Auto - Intense Bands)
Flat Field	Applied (Lens)
Serial Number	721BR00019
Software Version	2.0
Application	Colorimetric
Excitation Source	White epi illumination
Emission Filter	Standard Filter
Image Information	
Acquisition Date	9/18/2009 11:35:49 AM
User Name	Researcher
Image Area (mm)	X: 95.0 Y: 71.0
Image Pixels	X: 1392 Y: 1040
Pixel Size (um)	X: 68.2 Y: 68.2
Data Range (Int)	13665 - 64753
<div>OK</div> <div>Cancel</div>	

### ANALYSIS SETTINGS

The settings that were used when the gel was analyzed (for example, Band Detection and Molecular Weight Analysis, if they were performed) are displayed here.

### NOTES

You can add notes, make custom labels for each of the lanes, point out the type of samples used, or anything else you want to point out about the results.



## Displaying Data

The results from analyzed data associated with the gel images can be visualized using an Analysis Table, a Lane Profile, a Standard Curve, or as a Report.



Buttons to toggle these views on and off are located in the main toolbar. All of them can be viewed at one time. Each option is described below.

### Lane and band definitions

- Band Number—Each band in a lane has a unique number, sorted from top to bottom
- Band Label—User can assign a custom label to each band, by clicking in the Band Label field of the Lane and Band table
- Molecular Weight—The molecular weight of the band is calculated based upon the user-defined standard and regression method. Italic values indicate extrapolated values. When using nucleic acid gels, the size of the band is displayed in base pairs
- Relative Front—Values between 0-1 indicate the relative movement of the band from top to bottom
- Abs. Quant.—Absolute quantification of the band
- Rel. Quant.—Relative quantification of the band, compared to the reference band
- Band %—Percentage of the band volume compared to all band volumes in the lane
- Lane %—Percentage of the band's volume compared to the entire volume of the lane


### Volume definitions

- Volume Number—Each volume has a unique number assigned
- Volume Label—Software-generated labels for different types of volumes (U-Unknown, B-Background, S-Standard) can be changed by user in Volume Properties
- Volume—The sum of all the intensities within the band boundaries.
- Adjusted Volume —The background-adjusted volume
- Mean Background—The mean value of the background
- Absolute Quantity Volume—The quantity of the volume based on the standard volumes and the regression method
- Relative Quantity Volume—The ratio of the adjusted volume and the adjusted volume of the reference volume



- # Pixels—Number of pixels inside the volume boundary
- Minimum Value—Intensity of the pixel with the minimal intensity inside the volume
- Maximum Value—Intensity of the pixel with the maximum intensity inside the volume
- Mean Value—Mean value of all pixels inside the volume boundary
- Standard Deviation—Standard deviation of all pixels inside the volume boundary
- Area—Area of the volume in mm<sup>2</sup>

## Analysis Table Options

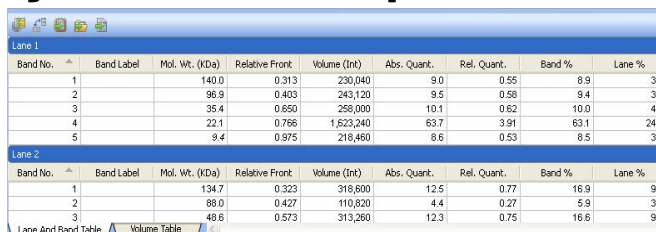

Band No.	Band Label	Mol. Wt. (KDa)	Relative Front	Volume (Int)	Abs. Quant.	Rel. Quant.	Band %	Lane %
1		140.0	0.313	230,040	9.0	0.55	8.9	3.5
2		96.9	0.403	243,120	9.5	0.58	9.4	3.7
3		35.4	0.650	258,000	10.1	0.62	10.0	4.0
4		22.1	0.766	1,623,240	63.7	3.91	63.1	24.9
5		9.4	0.975	216,460	8.6	0.53	8.5	3.3

Band No.	Band Label	Mol. Wt. (KDa)	Relative Front	Volume (Int)	Abs. Quant.	Rel. Quant.	Band %	Lane %
1		134.7	0.323	318,800	12.5	0.77	16.9	9.9
2		88.0	0.427	110,820	4.4	0.27	5.9	3.4
3		48.6	0.573	313,260	12.3	0.75	16.6	9.7

Numerical data associated with your analysis can be viewed in the analysis table. Data from the Lane and Band analysis can be viewed in the Lane and Band tab. If volume analysis was performed, this data can be viewed in the Volume Table tab.

Clicking on the Analysis Table button opens a tabular display of the data extracted from the results. The icons above the table offer several ways of displaying and exporting Analysis Table data.

## SET WINDOW SIZE

To change the size of your Analysis Table window, move your mouse to the top of the window until the cursor changes into a double-headed arrow. Left-click and drag the edge of the window until you can see all of your data.

**Note:** Resizing the Analysis Table window is restricted when a Protocol window is open.

## DISPLAY DATA OPTIONS

There are three tabs in this dialog box: Measurements, Display and Export.



**Measurements**—Choose the measurements you want displayed by clicking the arrow buttons to move them from the Displayed to the Not Displayed field.

**Default display settings**—“Move selected lane to top” is on or off, depending on whether the checkbox is selected or not.

**Per Measurement Precision**—You can set the precision (decimal places) for the measurements in both the Lane and Band table and the Volume table.

**Example Precision**—This field shows how your measurement will display, with the number of decimal places you have chosen in Per Measurement Precision.

**Export formatting**—This field offers two checkboxes to include lane headers (Lane and Band table tab only) and/or column headers in your exported file.

**Export delimiter**—This field offers three delimiter options for the exported file.

- Comma delimited CSV (Comma separated values)
- Tab delimited
- Other user-defined delimiter

## CHANGE ANALYSIS TABLE ORIENTATION

This button toggles between two table orientations.



**Horizontal**—Displays the lanes/volumes beside each other, so the user can scroll through the table from left to right.

**Vertical**—Displays the lanes/volumes on top of each other, so the user can scroll through the table from top to bottom.

## EXPORT AN ANALYSIS TABLE

The next three buttons offer several ways of exporting the data in the Analysis Table. See Chapter 10 for complete information on exporting your data.

**Copy Analysis Table to the Clipboard**—Click this button to copy the Analysis Table to your clipboard, in order to paste your Analysis Table to word processing or presentation applications. It is best to use the Vertical table orientation when copying to an 8-1/2x11” page, to allow the columns enough room to display.



**Export Analysis Table to a File**—When exported as a CSV (comma-separated values) file, your data file can be opened in a database application.




**Export Analysis Table to a Spreadsheet**—This function allows you to use Excel’s sorting and formula functions to manipulate your data. If you have this program installed on your computer, Excel (or Numbers, on a Mac) will open with your spreadsheet displayed.





## Lane Profile

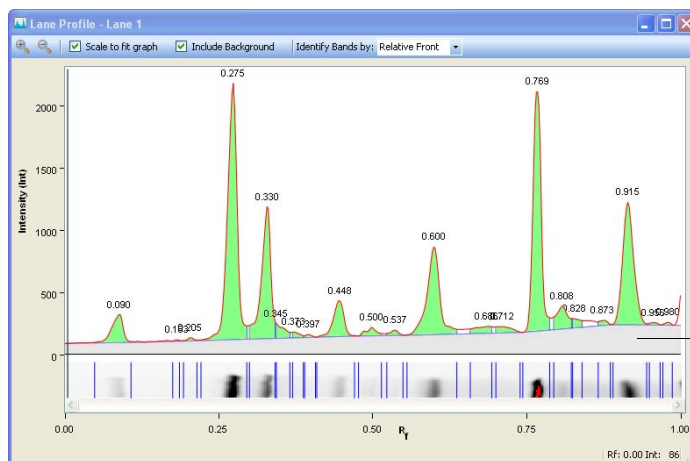

Background

The Lane Profile option shows you a cross-section of the selected lane, rotated 90 degrees. You can select other lanes by clicking on them while the Lane Profile window is open.

If the “Include Background” checkbox is selected, the Lane Profile window shows the subtracted background under the blue line, and the area used for band quantification in green under the red line.

You can see the current Rf (relative front) value and the intensity at the mouse cursor in the lower right corner of the Lane Profile screen.

You can change how the bands are labeled by choosing from the options in the dropdown list. By default, the bands are labeled with the band number.

### SCALE TO FIT GRAPH

You can choose the highest point of the display to define the range of the graph, to provide the best view of the lane profile.

You may want to clear the Scale to Fit Graph checkbox, to display the entire range of possible intensity values in your graph. This allows valid comparisons between different lanes.

### INCLUDE BACKGROUND

If you clear the Include Background checkbox, the area of the Lane Profile which represents the background of the image will not be displayed.





## IDENTIFY BANDS BY:

You can display callouts on the Lane Profile which indicate these attributes:

- Band Number
- Band Label
- Molecular Weight
- Relative Front
- Volume
- Abs. Quant.
- Rel. Quant.
- Band %
- Lane %

## Standard Curve

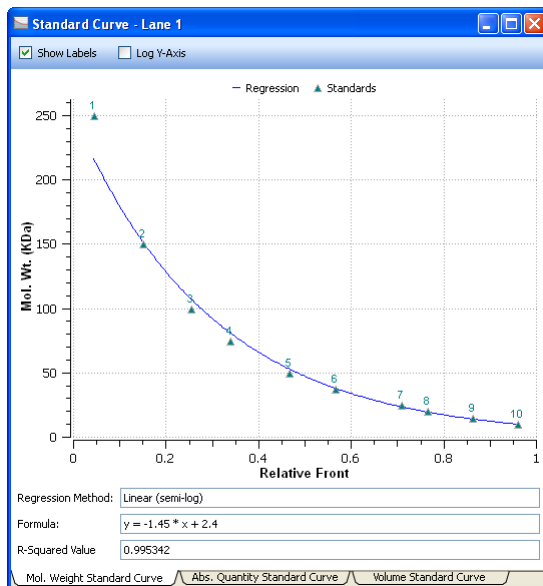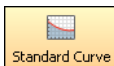

The Standard Curve dialog box displays the best curve fit for your defined standards and the bands relative to this curve, for the lane you have selected in the image. The tabs display the standard curves for three different analyses.



Standards are displayed in green, and unknown bands are in red. You can change the molecular weight display on the Y-axis between linear and log scale by clicking the Log Y-axis box at the upper left. The regression method you have chosen in Molecular Weight Analysis Tools is displayed, as well as the formula (if applicable) and the  $R^2$  value of the regression method.

Tabs in this window allow you to view the Mol. Weight Standard Curve, the Abs. Quantity standard Curve, or a Volume Standard Curve.



See Chapter 9 “Generating Reports” on page 79 to learn about the choices you have when deciding what to include in your report.








## 8 Analyzing Images

---

Analysis Tool Box tools are enabled once an image file is opened and in focus. An active or “in focus” window has a darker blue menu bar on a Windows PC. On a Mac, the window control icons display more brightly when a window is active. This distinction allows you to edit one of many open image files on your desktop.

### Analysis Tool Box Tools

#### Auto Analysis

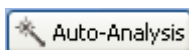


Clicking the Auto Analysis button in the Analysis Tool Box allows you to do the following:

- Analyze images that were obtained with protocols that did not include steps for detection and analysis
- Change your analysis parameters to re-analyze your images

**Note:** If you change any settings to an analyzed gel, the initial analysis will be overwritten. To preserve both analyses, save your image file with a different name.



## DETECTION SETTINGS

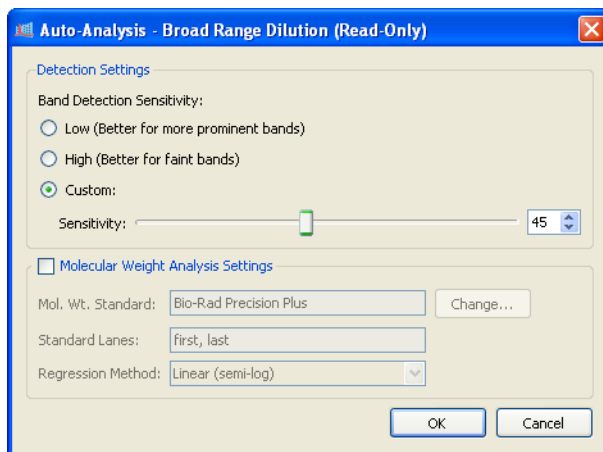

You have the following choices for band detection:

**Low Band Detection Sensitivity**— This setting will set detection at a low level, for images with more prominent bands. Faint bands will not be detected with this setting.

**High Band Detection Sensitivity**— This setting will set detection at a higher level, for images that are more faint. Extraneous bands can be removed later, using the Band Tools in the Analysis Tool Box. See “2. Lane and Band Tools” on page 62.

**Custom**— You can set a numerical value between 1-100, in order to select the best detection sensitivity for your sample.

**Note:** When Low or High Band Detection Sensitivity is used, these are the numerical values that are set:

Low sensitivity = 25

High sensitivity = 75



## MOLECULAR WEIGHT ANALYSIS SETTINGS

**Molecular Weight Standard**—Run any of the many Bio-Rad standards, or third-party standards which you have placed in your standards list. For information about adding standards to the standards list, refer to “Standards” on page 34. See “Ordering Information” on page 99 for catalog numbers.

**Standard Lanes**—Choose or change the lanes in which the standards are placed.

**Regression Method**—You have a choice of four regression methods. For complete information, see “Regression Methods” on page 36.

## Analysis Toolbox

All of the following tools customize the *analyzed* data in your image files. These tools are available only when an image file is open. Click a specific image to select among many windows which may be open on your desktop. The menu bar of the selected window becomes a darker blue than other windows.

Click on any of the Tool Box icons to access tools. Click the green arrow button to the right of the tool name to return to the Analysis Tool Box menu.


**Note:** Some of these tools delete the existing analysis.





## 1. Image Tools

To show the image view menu, click on the Image Tools button.

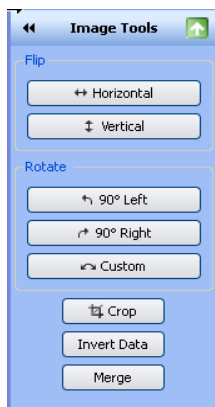

Click to return  
to the  
Analysis Tool Box  
menu

### FLIP

You can flip your gel image horizontally or vertically.

### ROTATE

You can rotate your gel image 90°, using the Left or Right buttons. To correct a slanted gel, click the Custom button and do the following:

1. Rotate the red arrows which appear over the gel to any degree between 0-360° by dragging them.
2. Right-click the gel image and choose Rotate, to set your gel in the new position. You will be prompted to Cancel or Continue, in case you want to reset the rotation.

### CROP

If you want only a portion of the gel image, follow these instructions to crop the image:

1. Click the Crop button.
2. Using the square handles, drag the red box to include the area you want to keep.
3. Right-click within the red box to remove the rest of the image. You will be prompted to Cancel or Continue, in case you want to reset the area.



## INVERT DATA

Used for negative stains and zymograms. Intensity values of bands must be greater than background, in order to perform analysis on the gel. View your gel as a 3-D projection to determine if the data needs to be inverted.

## MERGE

Use this button to merge a chemiluminescent blot image with a colorimetric image of the same blot. If you have used colorimetric pre-stained standards for your chemiluminescent blot, you can acquire an epi-white light image of the blot to show the standards, and a chemiluminescent image to show your immunodetection. These two images can then be merged into a combined image with both signals.

**Note:** Merging images can have an adverse effect on quantitation. If accurate quantitation is required, perform analysis on the original separate images. Only images of the same size can be merged.





## 2. Lane and Band Tools

Click the Lane and Band Tools button to select a tab, one to use the lane tools and one for band tools.

### LANE TAB

You can choose how your lanes are detected, using the Automatic or Manual Lane Finder buttons.



- Click the Automatic button if your gel image is fairly typical
- Click the Manual button if you only want to detect a specific number of lanes, or if automatic lane detection did not find all the lanes. You can then drag the handles on the corners of the red square, to resize the lane frame.



## ALL LANES

**Resize**—You can resize all of the lanes by dragging the handles on the corners of the red square to fit your gel image.

**Adjust**—If your gel image is somewhat irregular, you can adjust the orientation of all lanes by dragging a single corner of the lane frame.

You can also add additional anchor points on the top or bottom borders of the rectangle by clicking on the lane frame. Remove any unneeded anchor point by right-clicking it with your mouse. By dragging these anchor points, you can adjust for ‘smiling’ gels.

**Delete**—You can delete all lanes.

## SINGLE LANE

**Add**—You can add a single lane to your gel image. First click the Add button, then click within the lane frame where you want the new lane. The lanes will be renumbered.

**Note:** To add a lane outside the frame, add a lane inside the frame and use the move tool to expand the lane out of the frame’s boundaries.

**Delete**—You can delete a single lane. First click the Delete button, then click either on the lane or its lane number. The lanes will be renumbered.

**Bend**—You can bend a single lane to better fit your gel image. First click the Bend button, then click and drag one of the square anchors to fit your image.

You can add additional anchor points within the lane by left-clicking your mouse. Drag these anchor points to adjust the lane to fit the gel image. Remove anchor points by right-clicking them with your mouse.

**Move**—You can move a single lane to a new position on your gel image. First click the Move button, then click on the lane you want to move. Drag it to a new location. The lanes will be renumbered, according to their new position.

**Width**—You can change the width of a single lane.

## LANE BACKGROUND SUBTRACTION

Perform lane-based background subtraction by selecting Enable Subtraction in the Background Subtraction field. Use the Lane Profile view to see the subtracted lane background.

**Rolling Disk**—You can specify the size of a hypothetical “rolling disk” (between 1-99 mm) which removes background levels along the length of the lane. The size of the disk determines how much background will be subtracted.




A large disk will follow the profile trace less closely, touching fewer points along the trace and removing less background. A smaller disk will more closely follow the profile trace, removing more background.

A disk radius that is too large will result in poor background removal. A disk radius that is too small may subtract actual data. For most samples, a size of 10 mm or less is usually appropriate. You can perform this task several times until you are satisfied with the background removal.

**Apply to selected Lane**—If you select this box, the level of background subtraction specified with the above buttons will be applied only to the selected lane. This option gives you the opportunity to set different background subtraction levels for each lane.

## BANDS TAB

This tab allows you to detect bands or reset band detection settings. Click a button from the menu, then click in a lane of your image.

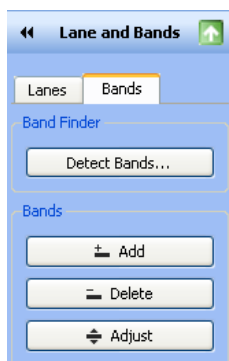

**Add**—Use this option to add faint bands. First click the Add button, then add a new band by clicking anywhere inside that lane.

**Note:** You can darken your entire image to view faint bands more easily by using the Gamma slider in the Image Transform dialog box. For complete instructions, see “Image Transform” on page 46.

**Delete** — You can delete bands that are not relevant to your analysis. First click the Delete button, and then click any bands you want to remove.



**Adjust**— You can adjust the height of a band. First click the Adjust button. Two boundary lines appear around each band. Move your mouse over a boundary line until you see a double-headed arrow. Move the boundary line up or down, the center is recalculated, and the band appears there.

**Note:** You can also adjust the band boundaries in the Lane Profile view.





### 3. Molecular Weight Analysis Tools

These tools enable you to determine molecular weight (or base pairs, if using nucleic acid gels) by comparing your test sample with known standards.

**Note:** You can view each band's molecular weight in the molecular weight column of the Lane and Band tab, in the Analysis Table view.

**MW Analysis Tools**

Standards  
Bio-Rad Precision Plus  
Change...

Standard Lanes  
first, last  
Select standard lanes by checking the box below the lanes

Regression  
Method:  
Linear (semi-log)

**Manage Standards**

Name	Type
Bio-Rad 20 bp Molecular Ruler	DNA
Bio-Rad 100 bp Molecular Ruler	DNA
Bio-Rad 100 bp PCR Molecular Ruler	DNA
Bio-Rad 500 bp Molecular Ruler	DNA
Bio-Rad 1 kb Molecular Ruler	DNA
Bio-Rad 2.5 kb Molecular Ruler	DNA
Bio-Rad AmpliSize Molecular Ruler	DNA
Bio-Rad Precision Molecular Mass Ruler	DNA

New  
Edit  
Copy  
Delete  
Rename

Show: DNA

Standards in red are built-in defaults.  
New standards can be created by copying a default.

OK



## MOLECULAR WEIGHT STANDARD

You can change the standards used to determine the molecular weight of your test samples. Click the Change button to access the Manage Standards dialog box, where you can choose another standard or add third-party standards.

## STANDARD LANES

Standard samples are placed in the first and last lanes, by default. You can specify other standard lanes by selecting the box below each lane, or by entering the standard lane numbers separated by commas. The Standard lanes are shown with the label “Std” below the lane.

## REGRESSION METHODS

There are four regression methods. See “Regression Methods” on page 36.




## 4. Quantity Tools

You can quantitate bands in your test samples automatically by using either the relative or absolute tabs under Quantity Tools.

## RELATIVE QUANTITY TAB

To compare the relative quantities of bands, you must first select the Quantity Tools Relative tab, as shown below.

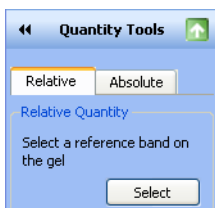


**To select a reference band from your image and quantitate all other bands, relative to the reference band:**

1. Click on the Select button.
2. Click on any band you want to use as a reference. A small “R” for Reference appears near the band you selected (see the red circle in the



upper left).

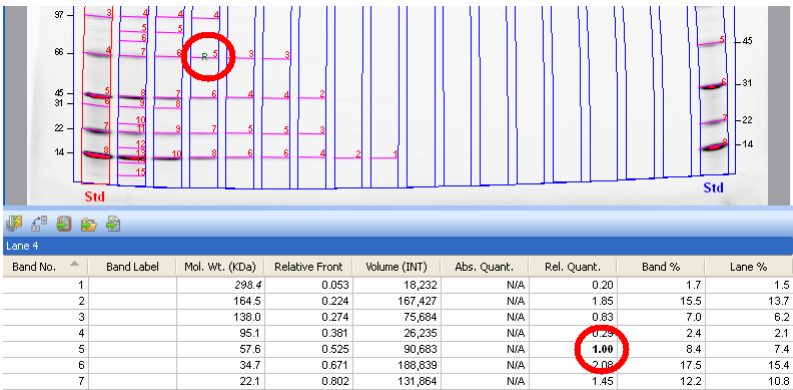

3. To review the relative band quantities, go to the Rel. Quant. column of the Analysis table, Lane and Band tab. The relative quantity is the ratio of the band volume divided by the reference band volume.

All of the other bands now display numerical values that are relative to your reference band. Values above 1.00 indicate that the band is more intense than the reference band. Values below 1.00 indicate the band is more faint.



## ABSOLUTE QUANTITY TAB

Absolute quantification is used to quantify bands based on known standard bands, using a calibration curve. To determine the absolute quantities of bands, first select the Quantity Tools Absolute tab, as shown below:

Quantity Tools

Relative Absolute

Absolute Quantity

Select Bands to add to calibration curve

Select

Standard Bands:

Ln	Bnd	Quantity
2	2	4.0
4	4	1.0

Delete

Units:

microgram

Regression

Method:

Linear

☐ Force through origin

Click on band to select a standard band

### To calculate the absolute quantities of the bands:

1. Click the Select button.
2. Select at least two standard (known) bands and assign quantity values. The values display in the Standard bands table. The greater the number of known bands, and wider the range of their values, the more accurate the absolute quantity calculation of the unknown bands will be.

**Note:** Any of these selections can be deleted. Select the entry in the Standard Bands field by clicking on it, and then click the Delete button below.



3. Select your unit of measure from the ten choices listed in the Units dropdown menu.
4. Select your regression method from the three listed in the dropdown menu. Keep the following guidelines in mind while making your choice.

**Linear**—This regression method generates a straight line that is the “best fit” of the values you provided, and is preferred in most cases.

**Point-to-Point**—This method generates a curve in which each data point is connected directly to the next, regardless of the shape of the resulting curve.

**Cubic spline**—This method generates a smooth curve which connects each data point. At least four standard points are required to use this method of least-squares polynomial fits.


Regression Method	Minimum number of standard bands	Minimum number with Force Through Option
Linear	2	1
Point-to-Point	2	1
Cubic Spline	5	4

5. Click on the Standard Curve button in the Toolbar, select the Abs. Quantity Standard Curve tab, and you will see a calibration curve. All of



your standards are represented by green triangles, and your Unknown values are represented by red triangles.

**Note:** Selecting the Force Through Origin checkbox will always start the standard curve graph at 0,0, regardless of the best curve fit.

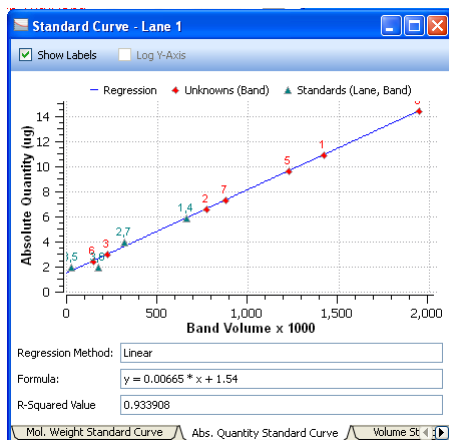


**Note:** Clicking any triangle in your Standard Curve table displays the numerical values associated with that band.





## 5. Annotation Tools

You can annotate your results with text and arrows, to draw attention to areas of interest.

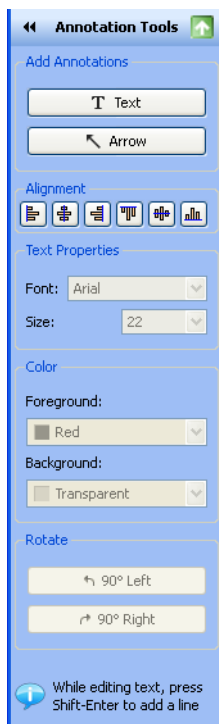

### ADD ANNOTATIONS

**Text**—You can add text annotations to your gel images to draw attention to important details. First click on the Text button, then on an area you want to emphasize. A text box appears with a dotted-line border. Type your comment into the text box. Click and drag the box to change its position.

**Note:** While editing text, press Shift-Enter to add text. This adds a new line in the annotation.

**Arrow**—To add an arrow, first click on the Arrow button. Click the area where you want the arrow to start, then drag to stretch the arrow point to the location you want to emphasize. To move the arrow, click on the middle and drag. To change where the arrow points, click on either end of the arrow. Square boxes appear; drag a box to change the length or orientation of the arrow.



## ALIGNMENT

**Alignment Icons**—These icons allow you to align multiple annotations, such as lane numbers, which you have manually added.

**To select annotations**—You can select several annotations by holding down the CTRL key and clicking each one, or by dragging a selection box around them.

**To copy annotations**—You can copy annotations within or between images. First select the annotation you want to copy. Then hold down the Ctrl key and press C to copy, and V to paste the annotations.

## TEXT PROPERTIES

You can change the size and type font of your text annotations. To select multiple items for change, hold down the keyboard Control button and left-click each item with your mouse.

**Font**—First click the text box you want to change. Then click the dropdown Font menu to show all fonts installed in your system. Click one of the font names to change the style of your text annotation.

**Size**—First click the text box you want to resize. Click the up and down Size arrows to increase or decrease the size of your text. You can set the font size between 6–72 points with the drop-down menu.

## COLOR

You can change the color of your text annotations to make them visible with any color scheme, and emphasize them further by adding a color to the annotation's background, which is invisible by default.

To change the color of multiple items, hold down the CTRL key and left-click each item.

**Foreground**—Click your text annotation or arrow. This enables the Foreground field, so you can use the drop-down menu to choose a different color.

**Background**—Click your text annotation. This also enables the Background field, so you can use the drop-down menu to choose a background color.

## ROTATE


You can rotate your text annotations 90 degrees to the left or right, by clicking the Rotate buttons.






## 6. Volume Tools

These tools give you a manual method for quantitating features on your sample image when automated lane and band analysis is not appropriate or possible, such as dot blots.

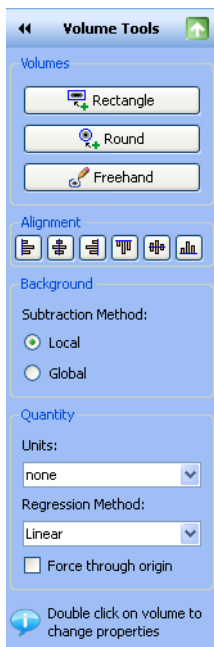

You can use the Volume Tools to quantitate the signal intensity of bands, spots, arrays, and other image data. You define an area of interest by surrounding it with a shape. You can choose a rectangle, a circle, or freehand shape by clicking the buttons under the Volumes field.

A default label appears within the shape you have drawn. The volume label can be one of three types; U for Unknown, Std for Standard, or B for Background, along with a number that is assigned sequentially.

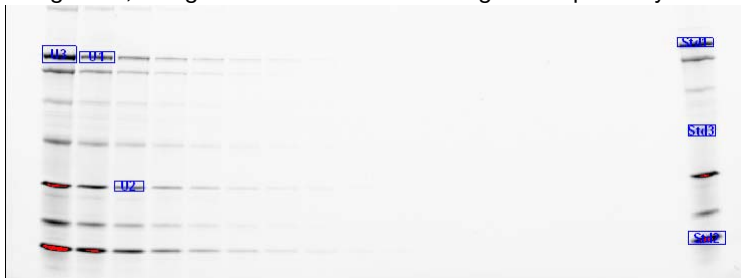



Each new volume you create initially has a red border, which indicates that the volume is selected. If you click elsewhere on the image, the border will change to blue, indicating that the volume is deselected

**Note:** Double-click a volume area to change its properties. You can define the volume type (Unknown, Standard, or Background), its quantity, or enter a custom name to replace the default label.



## RELATIVE VOLUME QUANTITY


You can choose any one volume as a Reference Volume by selecting the Reference Volume checkbox in the Volume Properties dialog box. The Reference Volume is indicated by an asterisk on the volume label, such as U1\*.

The relative quantities are displayed in the Relative Quantity column in the Volume Table. The relative quantity is the ratio of the background-adjusted volume divided by the background-adjusted reference volume.

All of the other volumes now display numerical values that are relative to your reference volume. Values above 1.00 indicate that the volume is greater than the reference volume. Values below 1.00 indicate the volume is less.



If you click the Standard Curve icon from the toolbar, a chart displays with all of the Unknown and Standard quantities.

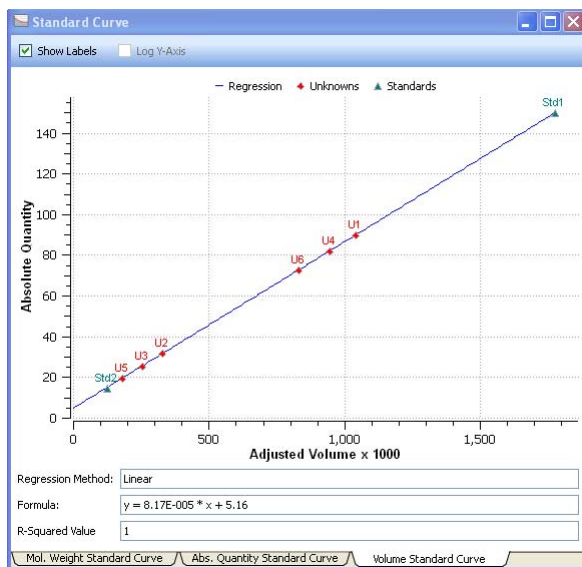

## ALIGNMENT

Align volumes by using the appropriate alignment icon. To select several volumes, CTRL-click each one or drag it to use a selection box.

Copy and paste selected volumes by pressing CTRL-C to copy and then CTRL-V to paste.

## VOLUME BACKGROUND SUBTRACTION

When you draw a volume, you will probably include some non-data background pixels inside the volume. These background pixels will usually have an intensity value that you do not want to include in your volume quantitation. There are two ways of calculating this background intensity for your volumes: local and global.



**Local**—Local background subtraction calculates a separate background intensity for each unknown and standard volume you create. For each volume, the intensities of the pixels in a 1-pixel border around the volume are added together and divided by the total number of border pixels. This gives an average intensity for the background around each, which is then subtracted from the intensity of each pixel inside the volume. If the background value is greater than the pixel value inside the volume, the background-adjusted quantity of the volume may be less than 0. In this case, you should redraw the border for this volume.

**Global**—Global background subtraction calculates a single background intensity for the entire gel. This average background intensity is then subtracted from all of the volumes in the gel.

### To calculate global background subtraction:

1. Create a volume using one of the volume tools in a representative background region of your image (that is, a non-data region similar to the background surrounding your data).
2. Double-click the volume. This opens the Volume Properties dialog box. Select the Background option button. The average intensity of the pixels in that background volume will be calculated and subtracted from each pixel in all standard and unknown volumes. Therefore, it is not necessary for the background volume area to be the same size as your unknown.

### Notes:

- If you select Global in the Volume toolbox but do not define a background volume as described above, no background subtraction will be performed
- If you create more than one background volume, all the pixels in those background volumes will be used to calculate the average background. Your background volume(s) will have default names B1, B2 . . . etc., based on the sequence they were created
- If the region you defined as background has a higher average intensity value than your data object, you will obtain a negative value for your adjusted volume in the Analysis Table. If this happens, select a new background region with less intensity than your data object

## VOLUME STANDARDS

If you have drawn your volume around an object of known quantity, you can use it to calculate the quantity of your unknown volumes.

1. To classify a particular volume as a standard, double-click it. This opens the Volume Properties dialog box.



2. Select the Standard option button, then enter the quantity in the quantity field. Click OK to close the dialog box. Your standard volumes will have default name S1, S2, etc., based on the sequence of their creation. The quantities of your unknown volumes is calculated based on the standard volumes and the selected regression method.

To review the regression curve, open the Standard Curve window and select the Volume Standard Curve tab.

## REGRESSION METHODS

Three regression methods are available to generate the volume quantitation curve: linear, point-to-point, and cubic spline. To display the standard curve, click on the standard curve icon in the tool bar and select the volume standard curve tab. See the Glossary to see how each of these methods is calculated.

The data for volume standards is found in the Absolute Quantity column of the Volume Table.

**Note:** Selecting the Force Through Origin checkbox will always start the standard curve graph at 0,0, regardless of the best curve fit.



## 9 Generating Reports


After viewing your results, you can generate a report that displays the analyzed gel images, all of the settings used in the protocol, and as much information about your data as you want to include.

The print report settings can be chosen with the Report Settings dialog box in the Edit menu, or from the Report button in the main toolbar.

### Report



Click the Report icon to produce a Print Preview of your report.

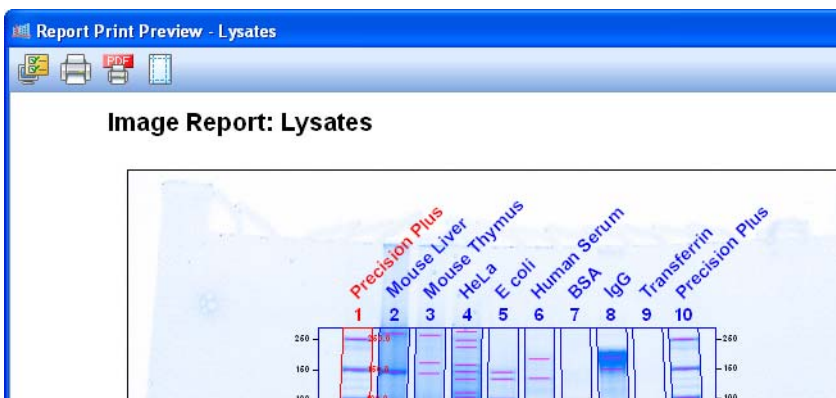





Clicking the Print icon prints all of the above information to a printer.


Use the following option screens to remove information from your reports. You are not deleting the data from your analysis, just information from your report.

## Display Report General Tab



The default setting is for all information to be included. These choices may have been made at other stages of your workflow, but you can customize your reports here based on the information you want to include.

Clear a checkbox to remove an item.

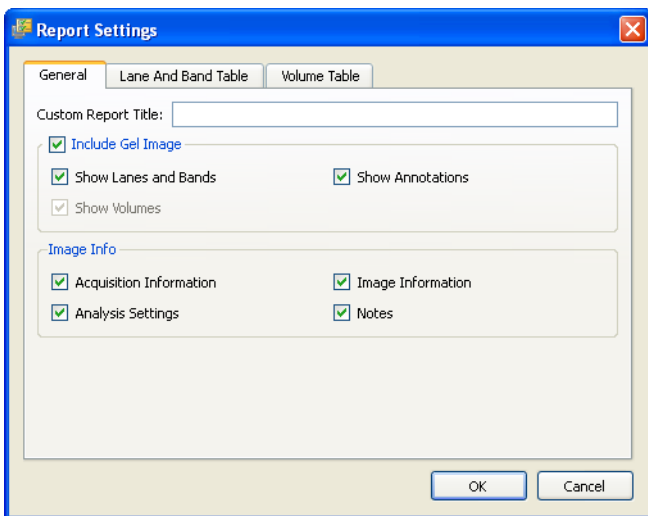

Clear any of these boxes to exclude the information from a report.

**Custom Report Title**—You can provide a custom report title for your report.


**Include Gel Image**—The following options will control which overlays display on your gel image.

- Show Lanes and Bands
- Show Volumes
- Show Annotations



**Image Info**

- Acquisition Information
- Analysis Settings
- Image Information
- Notes

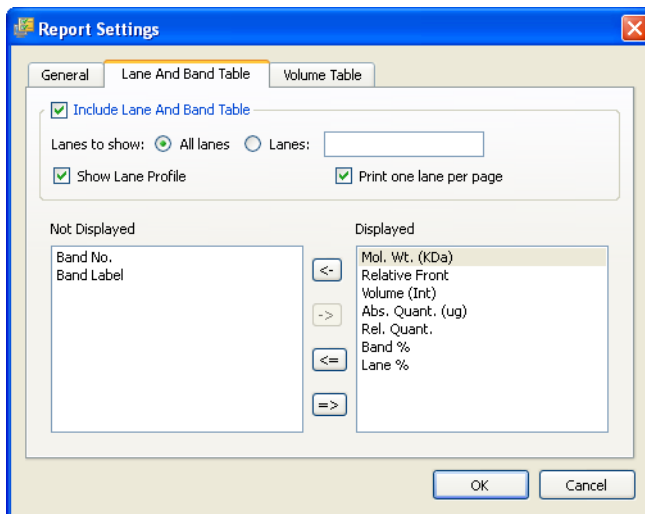**Lane and Band Table Tab**

Clear any of these boxes to exclude the information from your report.

**Include Lane and Band Table****Lanes to show:**


- All lanes or just user-defined lanes
- All lanes on one page, or one lane per page (add page break after each lane)

**Show Lane Profile**—Includes the lane profile for each lane.

**Move labels to the left** if you do not want their columns to be included in the report.



## Volume Table Tab

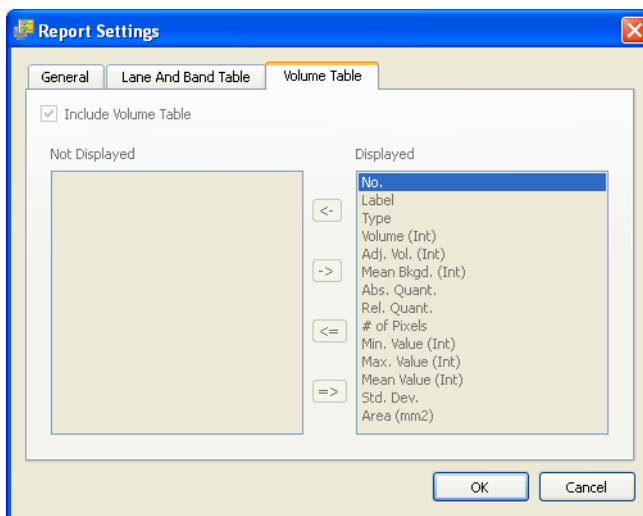

**Include Volume Table** - clear to exclude this information from your report.

**Move volume labels to the left** if you want to exclude their columns from the report.

## Print Report



Use the Print Report icon to print your report.

## Print Report to .pdf File



The Print Report to .pdf File icon brings up a Save dialog box so your .pdf file can be saved on your system.

## Adjust the Printer Settings



The Printer Settings icon provides options for paper size, orientation, and page margins.



# 10 Exporting Results

---

The most convenient way to archive complete information about your experiments is by producing reports, which is discussed in Chapter 9.

But you may want to export just your gel images or analysis table data for analysis in different programs, such as Quantity One®, FPQuest™, or InfoQuest™FP. Or you may need exported files for presentation or publication. This chapter describes all the ways Image Lab software can export your results.

## Exporting Gel Images


Image Lab software includes features for exporting gel images several ways.

- You can export displayed image data to a publication (choose Export for Publication)
- You can export raw image data as a 16-bit TIFF file (choose Export for Analysis)
- You can export image data to PulseNet; doing so reduces the image to an 8-bit TIFF file, limits its resolution, and restricts its file size to 300 Kb
- You can export lane and band tables as well as volume tables to a spreadsheet program or to a file

The options to export gel images are available from the Export option in the File Menu.

The available export options are described in the following sections.




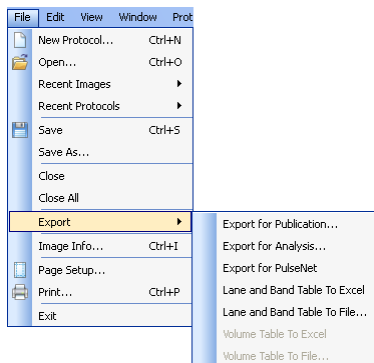

## Exporting Gel Images for Publication

Use this format only to export visual information to presentation or word processing software, such as PowerPoint or Word.


When you select File > Export > Export for Publication, you can export a displayed image to a file. You can select from .bmp, .png, .jpg, and TIFF formats. The gel displays with any lanes, bands, and annotations that appear on the screen.

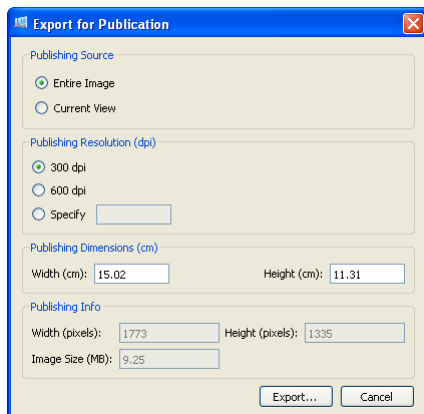


Note that you can select the entire image or the current view, select the resolution or specify a custom resolution, specify the publishing dimensions, and view the resulting published image size and dimensions.



You can zoom in on an area in a current view to export only that area or you can export the entire image. You can exclude annotations or overlays by

clicking the Display Gel Options  icon in the toolbar to access the appropriate settings.

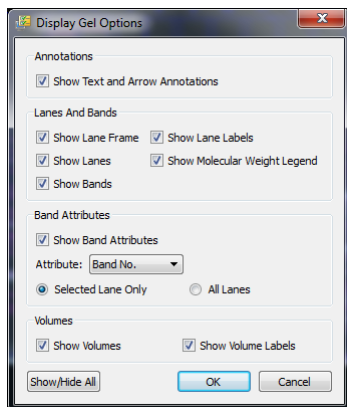

## Exporting Gel Images for Analysis

When you select File > Export > Export for Analysis, you can export full image data (raw data only) for analysis as a 16-bit TIFF file. The image may require contrast adjustment when it is imported into analysis software. This option creates a file that can be analyzed in other software such as Quantity One, FPQuest, or InfoQuestFP.

**Note:** Sixteen-bit TIFF images are not compatible with all image viewers.

## Exporting Gel Images to PulseNet

When you select File > Export > Export for PulseNet, Image Lab software reduces the image to an 8-bit TIFF image file. Resolution is limited and file size is restricted to 300 Kb.

## Exporting Lane and Band Tables to Excel

If you have Excel (or Numbers, on a Mac) installed on your computer, choosing File > Export > Lane and Band Table to Excel opens a table directly in Excel. You can then use the Save As option to produce other formats.



## Exporting Volume Tables to File

Choosing Export > Volume Table to File exports as a CSV file so the data file can be opened in a database application.

## Snapshot Tool Export

Use the Snapshot tool, available in the toolbar, to capture a displayed image to the clipboard or to save it to a file (.bmp, .gif, or .png).

## Analysis Table Export

You can export table analysis data from the File menu or by using the export buttons at the top of the Analysis Table window.

The Analysis Table window has several buttons to export data to different formats, depending on how the data are to be presented.

### COPY ANALYSIS TABLE TO THE CLIPBOARD



Click this button to copy the analysis table to your clipboard, and to paste your Analysis Table into word processing or presentation applications. It is best to use the Vertical table orientation when copying to an 8-1/2 x 11" page, to allow the columns enough room to display.

### EXPORT ANALYSIS TABLE TO A FILE



This exports as a CSV (comma-separated values) file, so your data file can be opened in a database application.

### EXPORT ANALYSIS TABLE TO A SPREADSHEET



This button allows you to use Excel's sort and formula functions with your data. If you have Excel (or Numbers, on a Mac) installed on your computer, your data opens in Excel.



# 11 Maintenance

---

This section includes instructions for maintaining the universal hood in proper working condition by replacing parts.

## UV Transilluminator Lamp and Starter Replacement

**Note:** The UV filter surface should always be kept clean from the chemical agents used as gel dyes. Use protective gloves when touching the UV transilluminator cover.

Depending on usage, the UV bulbs and starters last for many years. Replace bulbs when you notice them flickering. If a bulb does not turn on when it is new or moved, replace the bulb starter and test the bulb again.

There are three types of bulbs available. The catalog numbers are listed in Ordering Information [on page 95](#). The standard bulb is 302 nm. Optionally, the 254 nm bulb is used for cross-linking of protein, and the 365 nm is used to minimize denaturing of DNA.

### To replace the lamps:

1. Turn off the power.
2. Disconnect the power cord from the universal hood.
3. Remove the four screws located on the left/right sides of the transilluminator cover.
4. Remove the cover with the UV glass by sliding it forward, then lifting up.



5. Place it on a nonabrasive surface so that the glass does not get scratched or damaged.
- Note:** Do not put the UV cover directly on the bench. Wear gloves when touching the lamps.
6. Rotate the lamp until it becomes loose and the pins come to a vertical position.
  7. Remove the lamp. Install the new lamp by rotating so that the pins are horizontal and the lamp is tight.



8. Refer to the pictures above to replace a starter. Remove the starter by rotating the starter counterclockwise and then pull it out.
9. To replace a starter, insert it into the holder and rotate clockwise.
10. Reassemble the cover and retighten the screws on both sides.

## Epi-Illumination Lamp Replacement

The lamps are located behind two cover panels on the left and right sides of the universal hood's interior.






### To replace the lamps:

1. Turn system power off and remove the power cord.
2. Open the enclosure door.
3. Locate the epi-white light housing inside the hood and the screw on the center of the cover.
4. Use a socket wrench to remove the screw, and the plastic cover that it holds in place.
5. Pull the cover to remove it from the housing, and the lamp will become visible.
6. To remove the lamp, hold it at the plastic receptacle and then pull it from the receptacle.
7. Insert the new lamp into the lamp receptacle and then push it into the receptacle.
8. Reassemble the epi-white light cover.

## Fuse Replacement

Always unplug the instrument before changing or checking the fuses.

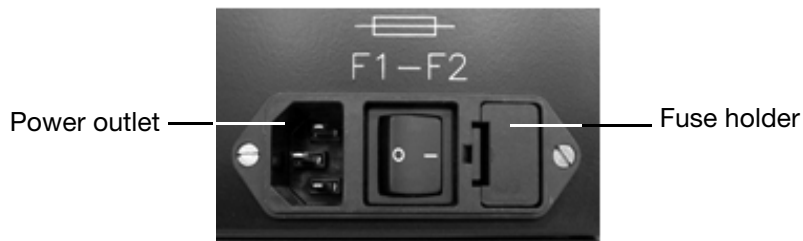

This unit is protected by 2 fuses (5 x 20 mm, 2A Slo-Blo). The fuses are located in the left side rear of the universal hood, in a fuse holder housed in the power entry module. See picture above.

### To replace the fuses:

1. Unplug the main power cable from the power outlet.
2. Use a screwdriver as a lever on the fuse holder to extract the fuse.
3. Remove the blown fuses and replace them with two new ones (catalog # 900-8935, quantity 10).
4. Slide the fuse holder into the power entry module until it snaps in place.







# Appendix A: Troubleshooting

Follow these suggestions to troubleshoot your Molecular Imager® system.

Problem	Possible Cause	Solution
Gel Doc XR+ and ChemiDoc XRS+ Systems		
Camera does not respond/camera not found	<ul style="list-style-type: none"> <li>• Camera power may be turned off</li> <li>• Camera cables may not be seated properly</li> <li>• Software driver for the camera is missing</li> <li>• Computer power-saving modes may be interfering with the camera driver</li> <li>• Cables may be defective</li> <li>• Camera may be defective</li> </ul>	<ul style="list-style-type: none"> <li>• Turn on the power to the camera</li> <li>• Make sure that all cables are connected as shown in the Installation Guide</li> <li>• If camera driver is not present, reload the camera driver from the Image Lab™ software CD</li> <li>• Disable power-saving modes on the computer</li> <li>• Replace cables</li> <li>• Replace camera</li> </ul>
Horizontal stripes in image when using the UV mode	<ul style="list-style-type: none"> <li>• Emission filter not in proper position</li> </ul>	<ul style="list-style-type: none"> <li>• Move filter lever so that the filter slider positions the filter under the camera lens</li> </ul>
Image is not visible on the monitor	<ul style="list-style-type: none"> <li>• Incorrect monitor settings</li> <li>• Lens cap attached</li> </ul>	<ul style="list-style-type: none"> <li>• See your computer manual for settings</li> <li>• Remove lens cap</li> </ul>
Image is not bright enough	<ul style="list-style-type: none"> <li>• Wrong emission filter in use</li> <li>• For chemiluminescent, emission filter is in front of lens</li> </ul>	<ul style="list-style-type: none"> <li>• Verify correct filter for application</li> <li>• Verify no emission filter is in front of lens</li> </ul>



Problem	Possible Cause	Solution
Printout does not look like the monitor image	<ul style="list-style-type: none"> <li>• Monitor settings are wrong</li> <li>• Printer settings are wrong</li> </ul>	<ul style="list-style-type: none"> <li>• Refer to monitor instructions to correct settings</li> <li>• Refer to the printer manual to correct settings</li> </ul>
Light leakage into the darkroom	<ul style="list-style-type: none"> <li>• The lens body is not seated properly against the gasket on the hood's adapter plate</li> </ul>	<ul style="list-style-type: none"> <li>• Loosen the thumbscrew and seat the lens properly against the gasket on the hood's adapter plate</li> </ul>
Unable to focus on the sample using white light transilluminator or conversion screen	<ul style="list-style-type: none"> <li>• Focus is not calibrated for samples using this light source</li> </ul>	<ul style="list-style-type: none"> <li>• Select Edit &gt; Instrument Setup to recalibrate focus using the proper accessory</li> </ul>
You are experiencing lens limits that seem artificially restricted	<ul style="list-style-type: none"> <li>• Camera lens not seated properly on lens mounting plate</li> </ul>	<ul style="list-style-type: none"> <li>• Reseat camera on lens mounting plate</li> </ul>
<b>Gel Doc XR+ System Only</b>		
Hot (very bright) pixels are seen in the image (Gel Doc XR+ system only)	<ul style="list-style-type: none"> <li>• Corrupted dark image correction</li> </ul>	<ul style="list-style-type: none"> <li>• In Edit &gt; Instrument Setup, recalibrate the dark image correction file</li> </ul>
<b>ChemiDoc XRS+ System Only</b>		
Camera does not respond/camera not found	<ul style="list-style-type: none"> <li>• Multiple FireWire devices are installed on the same computer (ChemiDoc XRS+ with FireWire only)</li> </ul>	<ul style="list-style-type: none"> <li>• Cycle power: first turn off the camera, hood and computer; then turn on camera, hood and computer, in that sequence</li> </ul>
ChemiDoc XRS+ images are grainy	<ul style="list-style-type: none"> <li>• Low signal-to-noise</li> </ul>	<ul style="list-style-type: none"> <li>• For chemiluminescent applications, make sure no filter is used</li> <li>• For applications other than chemiluminescent, make sure the filter is in line with the lens</li> <li>• Use sample that generates more signal</li> <li>• Use longer exposure times</li> </ul>



# **Appendix B: Accessories**

---

## **Calibrating Accessories**

If you are installing accessories along with your original system installation, you calibrate your system with a one-time Instrument Calibration wizard. Complete instructions are in the Installation Guide which arrives with your system, either document 10016995 for Gel Doc XR+ or document 10016999 for ChemiDoc XRS+. If you acquire new conversion screens, light sources or filters for a Gel Doc XR+ or ChemiDoc XRS+ imaging system after your original system installation, you will need to recalibrate your system to use them.

Refer to Chapter 4 for instructions on how to calibrate newly-acquired accessories.

## **Installing Optional Accessories**

### **UV/White Light Conversion Screen**

This optional white light conversion screen converts the UV light generated in the universal hood to white light. Your system must be calibrated to use the White Light Conversion Screen.

To calibrate, go to the Edit > Instrument Setup dialog box and select the appropriate checkbox. The Image Lab software will guide you through the calibration.



**To use the White Light Conversion Screen:**

1. Center the conversion screen on the imager stage.
2. Center your samples on top of the conversion screen.



3. Image the gel using your preferred application.


**XcitaBlue™ Conversion Screen**

The optional XcitaBlue™ conversion screen kit (catalog #170-8182) converts



UV to blue light, which allows you to visualize DNA samples while protecting against UV damage.

The XcitaBlue conversion screen is held in place by adhesive-backed edge guides. After installing the edge guides, the conversion screen remains centered and will not slide, even if you close the drawer rapidly.

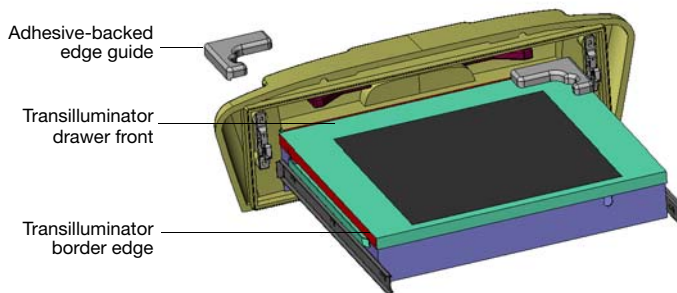




### To install the XcitaBlue Conversion Screen:

1. Perform a trial placement first, without removing the paper tape. Place the edge guides in each corner of your transilluminator, as shown above. The edge guides should touch the inside of the drawer front and fit over the edge of the metal transilluminator border (shown in red).
2. Remove the paper tape from the bottom surface of each edge guide.
3. Press each edge guide into position carefully, as it is difficult to reposition them once the adhesive surfaces touch.
4. Calibrate your system to use this accessory, by going to Edit > Instrument Setup. Select the “XcitaBlue Conversion Screen” checkbox under Illumination Options. The software will prompt you to calibrate the focus with height offset.
5. When you need to visualize a sample using the XcitaBlue conversion screen, place the screen between the edge guides.
6. Center the gel on top of the XcitaBlue conversion screen, and proceed with normal image capture. See the Gel Alignment Template Kit on the next page, for help centering your gels easily and consistently.

## Gel Alignment Template Kit

The Bio-Rad gel alignment template kit (catalog #170-8184) allows four sizes of standard agarose gels to be centered quickly and easily, as well as ensuring the consistent placement of each gel.

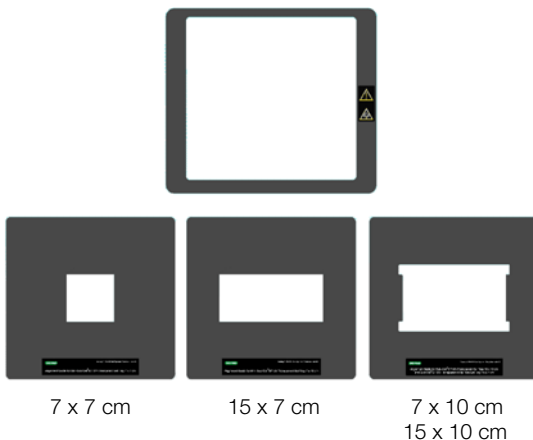

The kit contains:

- Magnetic locator frame



- Instruction sheet
- Alignment guides for the following gel trays:
  - Sub-Cell® GT UV-transparent mini-gel tray, 7 x 7 cm
  - Sub-Cell GT UV-transparent wide mini-gel tray, 15 x 7 cm
  - Sub-Cell GT UV-transparent mini-gel tray, 7 x 10 cm
  - Sub-Cell GT UV-transparent gel tray, 15 x 10 cm

The gel alignment templates can be used with the XcitaBlue™ conversion screen kit (catalog #170-8183). The templates fit exactly into the XcitaBlue conversion screen frame.

**To install and use the Gel Alignment Template Kit:**

1. Place the locator frame over the transilluminator, with the magnetic side down. Match the corners of the magnetic locator frame with the edges of the transilluminator. The UV symbol on the frame will be in the same orientation as the UV symbol on the imager.
2. Place the gel alignment template that matches the size of your sample tray or agarose gel into the magnetic locator frame.
3. Place your gel or gel tray into the open area of the template.

**Note:** No recalibration is necessary to use the Gel Alignment Template Kit.



## White Light Transilluminator

The white light transilluminator provides a white light that is independent of the UV bulbs. Between uses, this transilluminator can be stored in the rear of the universal hood.

### To install the white light transilluminator:

1. Open the enclosure door and pull the drawer toward you.
2. Locate the power cable positioned on the inside left side of the enclosure, behind the drawer slide.
3. Remove the black rubber boot that covers the connector plug and insert it into the white light transilluminator. Make sure that the main power switch on the transilluminator is in the ON position.

**Note:** Power to the white light transilluminator is controlled from either the software or the membrane touch pad on the universal hood.



4. Place the white light transilluminator on top of the UV transilluminator.



Your system must be calibrated to use the White Light Transilluminator.



To calibrate, go to the Edit > Instrument Setup dialog box and select the appropriate checkbox. The Image Lab software will guide you through the calibration.

## Orange Fluorescence Reference Plate

The Orange Fluorescence Reference Plate allows you to apply UV flat fielding corrections to your Gel Doc XR+ and ChemiDoc XRS+ imagers. Corrections are made for all UV illumination sources, filters, and the camera lens.

The Orange Fluorescence Reference Plate can be used to correct for image non-uniformities for red and orange gels, including:

- Ethidium bromide
- GelRed
- Flamingo
- Coomassie Fluor Orange
- SYPRO Ruby
- Krypton
- Qdots 625

### USE IT WITH:



The Orange Fluorescence Reference Plate (catalog #170-8008) can be used with several Bio-Rad Molecular Imager® systems, including:

- Gel Doc XR+ with Image Lab software
- ChemiDoc XRS+ with Image Lab software
- ChemiDoc XRS+ with Quantity One

When you order the reference plate you receive:

- Orange Fluorescence Reference Plate, overall dimensions 29.5 x 29.5 cm; viewing surface, 27 x 27 cm
- instruction sheet (10017296)



## CALIBRATING YOUR SYSTEM TO USE THE ORANGE FLUORESCENCE REFERENCE PLATE

Your system must be calibrated to use the Orange Fluorescence Reference Plate.

To calibrate, go to the Edit > Instrument Setup dialog box and select the appropriate checkbox. The Image Lab software will guide you through the calibration.

## Ordering Information

The following table contains catalog numbers and descriptions for all parts available for the Molecular Imager Gel Doc XR+™ or ChemiDoc XRS+™ systems, plus all optional accessories and replacement parts. For complete information, see the Bio-Rad catalog.

Catalog #	Description
<b>Molecular Imager System (includes Universal Hood, camera, cables, and accessories)</b>	
170-8195	Gel Doc XR+ with Image Lab System for PC/Mac
170-8265	ChemiDoc XRS+ with Image Lab System for PC/Mac
<b>Installation Kits</b>	
170-8199	Gel Doc XR+ Installation Kit
170-8299	ChemiDoc XRS+ Installation Kit
<b>Universal Hood</b>	
170-8180	Universal Hood
<b>Imaging Cameras</b>	
170-8181	Gel Doc XR+ camera with motorized zoom lens
170-8255	ChemiDoc XRS+ camera with motorized zoom lens
<b>Image Lab Software</b>	
170-9690	Image Lab software, Windows/Mac
<b>Optional Filters</b>	
170-8074	Filter, 520DF30, GFP
170-8075	Filter, 560DF50, Cy3/Rhodamine
170-8076	Filter, 630BP30, SYPRO Ruby
<b>Optional Accessories</b>	
170-8001	UV/White light conversion screen (UV to white light)
170-7950	White light transilluminator (Universal Hood plug-in)



<b>Catalog #</b>	<b>Description</b>
170-8182	XcitaBlue (UV to blue light) conversion screen kit, without standard detection filter
170-8183	XcitaBlue (UV to blue light) conversion screen kit, with standard detection filter
170-8008	Orange Fluorescence reference plate
170-8089	Printer, digital, Mitsubishi P93DW, 100/240 V, USB
170-3759	Bio-Rad fluorescent ruler
170-3760	Gel cutter ruler
170-8184	Gel Alignment Template Kit

### Replacement Parts

170-8026	Focus Calibration Target
170-8027	Image Lab Flat Fielding Disc
170-8185	XcitaBlue Viewing Goggles
170-7581	Mitsubishi thermal printer paper, 4 rolls
170-7813	Sample holders for gels
100-2784	UV transilluminator lid (includes UV filter glass)
1001-4106	Thumbscrew for camera
170-8081	Standard emission filter replacement, 62 nm
100-1370	UV bulb starter, quantity 3
930-2242	Multicolor target
<b>Lamps</b>	
100-2827	Lamp, epi-illumination, 5 W
100-1361	UVB Lamp, 302 nm (1 each)
170-8097	302 nm Lamp Kit, (6 lamps)
170-8098	254 nm Lamp Kit, (6 lamps)
170-6887	365 nm Lamp Kit, (6 lamps)
<b>Fuses</b>	
900-8935	Fuse T 2 A, 250 V, quantity 10
900-0234	Fuse T 4 A, 250 V, quantity 10
<b>Universal Hood</b>	
100-2785	Universal Hood right shield epi-illumination (lid complete)
100-1948	Universal Hood opal filter epi-white illumination (without metal frame)
100-2786	Universal Hood left shield epi-white illumination (lid complete)
100-2787	Universal Hood feet, 4
170-8068	Shield, UV, Universal Hood



Catalog #	Description
<b>Connection Cables</b>	
931-0071	Cable, USB, Type A to B, 10 ft.
901-0064	Cable, USB, Type A to B, 6 ft.
<b>Optional Analysis Software</b>	
170-9600	Quantity One Software
170-9300	FPQuest Software
170-9310	InfoQuestFP Software
<b>Protein Standards</b>	
161-0363	Precision Plus Unstained Standards
161-0373	Precision Plus All Blue Standards
161-0374	Precision Plus Dual Color
161-0375	Precision Plus Kaleidoscope
161-0385	Precision Plus WesternC Pack
161-0318	Prestained SDS-PAGE Standards, Broad Range
161-0317	Unstained SDS-PAGE Standards, Broad Range
<b>Nucleic Acid Standards</b>	
170-8351	EZ Load 20 base pairs Molecular Ruler
170-8352	EZ Load 100 base pairs Molecular Ruler
170-8353	EZ Load 100 base pairs PCR Molecular Ruler
170-8354	EZ Load 500 base pairs Molecular Ruler
170-8355	EZ Load 1 kb Molecular Ruler
170-8205	2.5 kb Molecular Ruler
170-8200	AmpliSize Molecular Ruler
170-8356	EZ Load Precision Molecular Mass Ruler (base pairs/ng of sample)
<b>Pulsed Field Standards and Markers</b>	
170-3624	CHEF DNA Size Standard, 5 kb ladder
170-3707	CHEF DNA Size Standard, 8-48 kb
170-3635	CHEF DNA Size Standard, lambda ladder
170-3605	CHEF DNA Size Marker, 0.2-2.2 Mb
170-3667	CHEF DNA Size Marker, 1-3.1 Mb
170-3633	CHEF DNA Size Marker, 3.5-5.7 Mb







# **Appendix C: Mitsubishi P93/P95 Thermal Printer Settings**

---


## **Windows**

The printer driver can be found on the Image Lab™ software installation CD in the Misc directory.

### **To set up a thermal printer on a Windows system:**

1. Install the printer driver.
2. Open the printer section in Control Panel.
3. Click the thermal printer icon and select Printing Preferences.



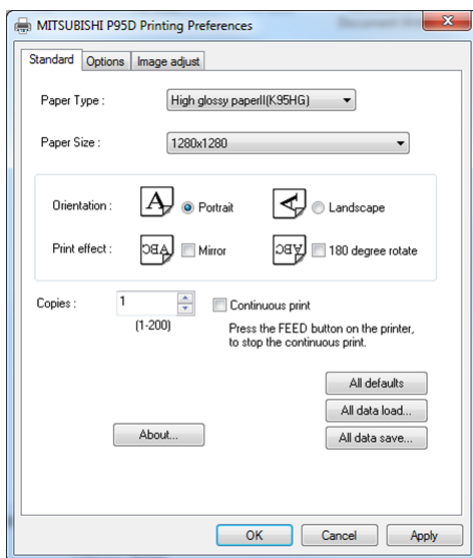

4. Configure the correct paper size: select 1280 x 1280.
5. Select OK to apply your changes.

## Macintosh

The printer driver can be found on the Image Lab™ software installation CD in the Misc directory.


### To set up a thermal printer on a Macintosh system:

1. Install the printer driver.
2. Connect the printer to the computer.

### To configure the correct paper size:

1. Start Image Lab software.
2. Select File > Page Setup.



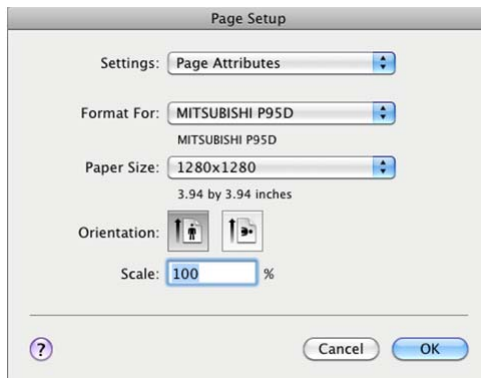

3. In the Settings drop-down menu, select Page Attributes.
4. In the Format For drop-down menu, select the Mitsubishi printer.
5. In the Paper Size drop-down menu, select 1280 x 1280.
6. In the Settings drop-down menu, select Save as Default.
7. Click OK to save the settings.







## Appendix D: Regression Calculation Methods

---

Each regression method calculates a standard curve. Some of the methods provide the formula for the standard curve. In this case, the molecular weight can be calculated by:

$x$  = relative front of the band of interest

$y$  = molecular weight of the band of interest

**Linear (semilog):** The linear equation is  $y = a + bx$ , where  $a$  is the intercept and  $b$  is the slope of the line.

**Note:** The linear equation is calculated on the **log** of the molecular weight values.

The  $R^2$  value may be used to determine the overall quality of the linear fit. A linear regression with an  $R^2$  value of  $>0.99$  is considered a very good fit. The primary advantage of this method is that it is extremely simple. The primary disadvantage is that it will deliver incorrect results if the data are not very linear.

**Point-to-Point (semilog):** No single equation is available for the point-to-point method. The slope of each segment of the curve between data points is calculated independently.

**Note:** The log of the molecular weight values is used to calculate the slope for each segment of the curve.



**Logistic:** The Logistic-4PL equation is  $y = d + \frac{a - d}{1 + \left(\frac{x}{c}\right)^b}$

where:

$x$  is the mobility

$y$  is the molecular weight

$a$  is the estimated molecular weight at infinity

$b$  is the slope of the tangent at midpoint

$c$  is the midpoint

$d$  is the estimated molecular weight at zero mobility

Since the curve generated by the logistic-4PL regression method represents a perfectly shaped S, it might not fit the data very well in all cases.

**Cubic Spline:** Cubic spline curves are smooth curves that go through every data point. The model is a cubic polynomial on each interval between data points. In some cases, a spline curve can work well as a standard curve for interpolation. However, because the curve is calculated individually for every pair of points, it does not correspond to any single equation.



# Glossary

---

**CCD:** (Charge-coupled device) A light-sensitive silicon chip used as a photodetector in Gel Doc XR+ and ChemiDoc XRS+ camera systems.

**Colormaps:** Different color representations of a gel image

**Electrophoresis:** A technique for separating molecules based on the differential movement of charged particles through a matrix when subjected to an electric field.

**Example precision:** The number of decimal places chosen for displaying a measurement.

**Flat Fielding:** An average intensity computation which compensates for non-uniformities generated by an instrument.

**Histogram:** A graphed representation of the brightness, or gray value, of an image.

**Native charge density:** The inherent electrical charge of a protein without the addition of SDS.

**pI:** Isoelectric point; the pH at which a protein molecule carries no net charge.

**Rf:** Relative front value of the band. In Image Lab software, Rf has a value between 0-1, and indicates the relative movement of the band from top to bottom.



**Quantitative Imaging:** Determines the quantity of a protein's components, through analysis of the pixel values in a digital image of the sample.

**UV-B:** The range of ultraviolet light used by the system.

**UV transilluminator:** The part of the imager that transmits UV light through a sample.



# Index

---

## A

Accessories 93  
Analysis Table  
    change orientation 52  
    copy table to the clipboard 52, 86  
    export to a file 52, 86  
    export to a spreadsheet 52, 86  
Analysis Toolbox  
    Annotation Tools 72  
    Auto Analysis 57  
    Image Tools 60  
    Lane Tools 62  
    Mol. Wt. Analysis Tools 66  
    Quantity Tools 67  
    Volume Tools 74  
Application Table 40

## B

Band detection 62

## C

Calibration  
    initial installation 15  
CCD  
    defined 109  
CCD camera and Universal Hood 2  
CCD Cameras 2  
ChemiDoc XRS+  
    product description 2  
    workflow 11  
ChemiDoc XRS+ Applications 13  
Choosing a Protocol 37  
Computer Requirements 4  
Computer Specifications 5  
Conversion screens 4  
Creating Custom Protocols 27  
Cubic spline  
    defined 108

## D

Display  
    Gel Options 44



Display Data Options 51

## **E**

Editing Protocols 38

Electrophoresis

defined 109

Emission filters 3

Epi-Illumination Lamp Replacement  
88

Exporting results 83

## **F**

Fit in Window

how to resize Results Files 45

Flat Fielding

defined 109

Fuse Replacement 89

Fuses iii

## **G**

Gel Doc XR+

product description 1

Gel Doc XR+ Applications 8

Gel Doc XR+ Imager Workflow 7

Generating Reports 79

## **I**

Image Colors

colormap choices 48

how to change colormaps 47

Image Lab 2.0 Software 3

Image Transform

how to set image contrast 46

## **L**

Lane detection 62

## **M**

Maintenance 87

## **O**

Optional Accessories

installing 93

Optional conversion screen 4

Optional Printer 4

Ordering Information 99

## **P**

Power Safety ii

Protocol File

compared to a Results File 17

Protocol Files defined 17

Protocols, creating custom 27

Protocols, Opening and Editing 38

## **R**

Report

adjusting printer settings 82

printing 82

printing to .pdf file 82

Reports 79

generating 79

Results

exporting 83

Results File

compared to Protocol File 17

Results Files



- overview 43
- viewing as a 3D projection 48
- viewing as a Lane Profile 53
- viewing as a Report 55
- viewing as a Standard Curve 54
- viewing as an Analysis Table 51
- viewing Image Information 49

Running a Protocol 39

## **S**

Safety Information i

System Calibration 15

## **T**

Technical Specifications

- ChemiDoc XRS+ 14
- Gel Doc XR+ 9

Three-dimensional projection 48

Troubleshooting 91

- ChemiDoc XRS+ only 92
- Gel Doc XR+ only 92

## **U**

Universal Hood

- product description 3

UV Starter replacement 87

UV Transilluminator replacement 87

## **V**

venience 4

Viewing Images 43

Voltage Setting ii

Volume Tools 74

## **W**

Warranty i

White Light conversion screen

- description 4

White Light Transilluminator 4

Workflow

- Step 1. Gel Imaging 28
- Step 2. Detect Lanes and Bands 33
- Step 3. Analyze Molecular Weight 33
- Step 4. Specify Content of Reports 37

## **X**

XcitaBlue Conversion Screen 4

## **Z**

Zoom tools

- how to resize images 45





**Bio-Rad  
Laboratories, Inc.**

Life Science  
Group

Web site [www.bio-rad.com](http://www.bio-rad.com) **USA** 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 01 877 89 01 **Belgium** 09 385 55 11 **Brazil** 55 31 3689 6600  
**Canada** 905 364 3435 **China** 86 20 8732 2339 **Czech Republic** 420 241 430 532 **Denmark** 44 52 10 00 **Finland** 09 804 22 00 **France** 01 47 95 69 65  
**Germany** 089 31 884 0 **Greece** 30 210 777 4396 **Hong Kong** 852 2789 3300 **Hungary** 36 1 459 6100 **India** 91 124 4029300 **Israel** 03 963 6050  
**Italy** 39 02 216091 **Japan** 03 6361 7000 **Korea** 82 2 3473 4460 **Mexico** 52 555 488 7670 **The Netherlands** 0318 540686 **New Zealand** 0508 805 500  
**Norway** 23 38 41 30 **Poland** 48 22 331 99 99 **Portugal** 351 21 472 7700 **Russia** 7 495 721 14 04 **Singapore** 65 6415 3188 **South Africa** 27 861 246 723  
**Spain** 34 91 590 5200 **Sweden** 08 555 12700 **Switzerland** 061 717 95 55 **Taiwan** 886 2 2578 7189 **United Kingdom** 020 8328 2000